




OPTILAB

OP301

Portable Optics Laboratory

Product Manual

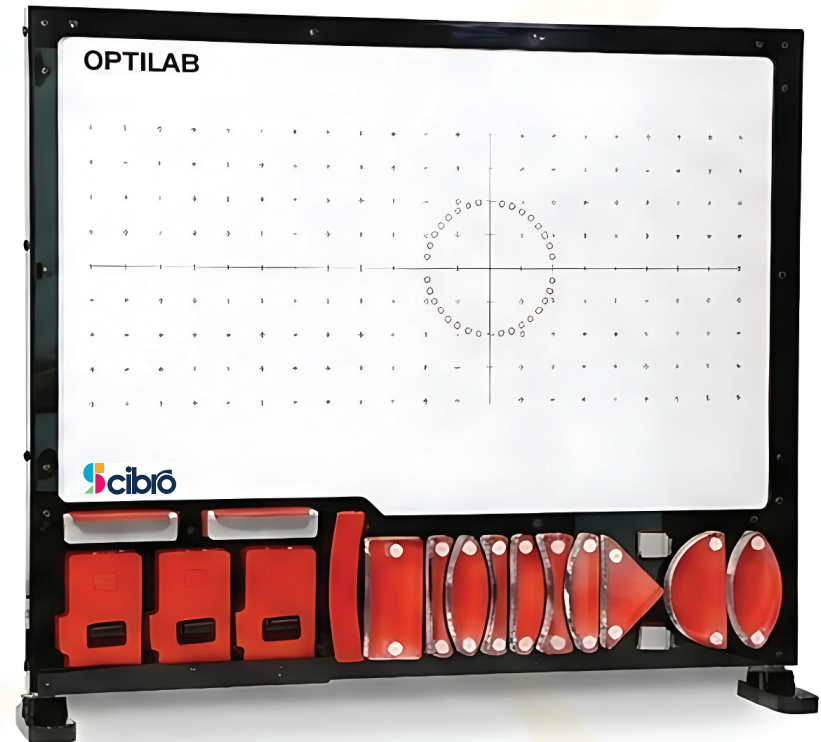

www.scibro.com

CONTENTS

LASER OPTICS SET OP301 – Portable Optics Laboratory

Part 1: Plane Mirror

Angle of incidence	4
Rotation of mirror and reflection angle	5
Perpendicular Mirrors	6
Parallel Mirrors	7
Periscope (Submarine Eye)	8
Virtual Image from Real Source	9
Virtual Image (Hypothetical light source behind the mirror)	10

Part 2: Concave Mirror

Mirror Vertex	11
Mirror Center	12
Principal Focal Point	13
Real Image (Magnification equal to unity).....	14
Real Image (Infinite rays)	15
Real Image (Infinite sources)	16

Part 3: Convex Mirror

Mirror Vertex	17
Mirror Center	18
Principal Focal Point	19
Virtual Image (Magnification less than unity).....	20

Part 4: Refraction of Light

Transmission between two media	21
Total reflection & critical angle	22

Part 5: Plane-Parallel Slab

Transmission through slab	23
---------------------------------	----

Part 6: Prism

Deflection Angle	24
Total Reflection (90°)	25
Two Total Reflections (180°)	26

Part 7: Convex Lens



Optical Center	27
Principal Focal Point	28
Effect of Convexity	29
Real Image (Magnification = 1)	30
Real Image (Infinite rays)	31

Part 8: Concave Lens

Optical Center	32
Virtual Focal Point	33
Virtual Image	34

Part 9: Physics of the Eye

Healthy Lens	35
Myopia (Correction)	36

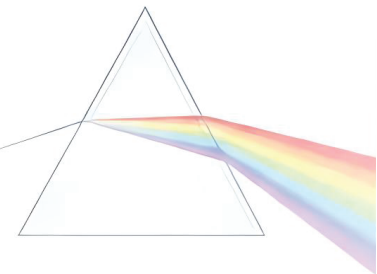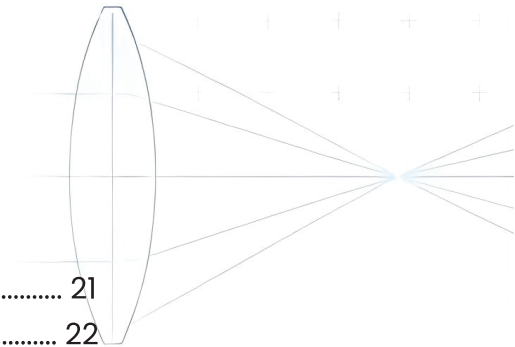


LASER OPTICS SET

OP301

Portable Optics laboratory.

Easy and interesting way to studying geometrical optics based on experiments.

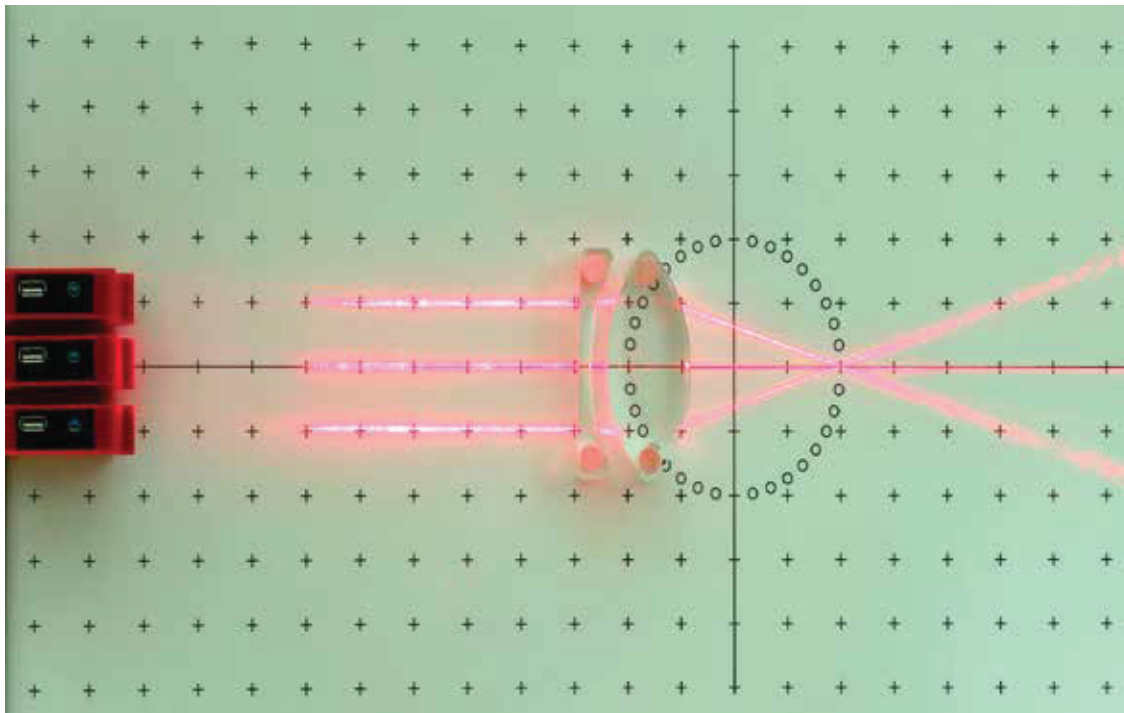


List of experiments (33 Experiments)

Part 1: Plane Mirror	(7 Experiments)
Part 2: Concave Mirror	(6 Experiments)
Part 3: Convex Mirror	(4 Experiments)
Part 4: Refraction of Light	(2 Experiments)
Part 5: Plane-Parallel Slab	(1 Experiments)
Part 6: Prism	(3 Experiments)
Part 7: Convex Lens	(5 Experiments)
Part 8: Concave Lens	(3 Experiments)
Part 9: Physics of Eye	(2 Experiments)

1- Plane Mirror

1-Angle of incidence


- Place the plane mirror on the determined position in the figure.
- Place the light source and illuminate the plane mirror along the **guiding line**.
- By varying the incidence angle, study the **relation between the incidence and reflection angles** in the plane mirror

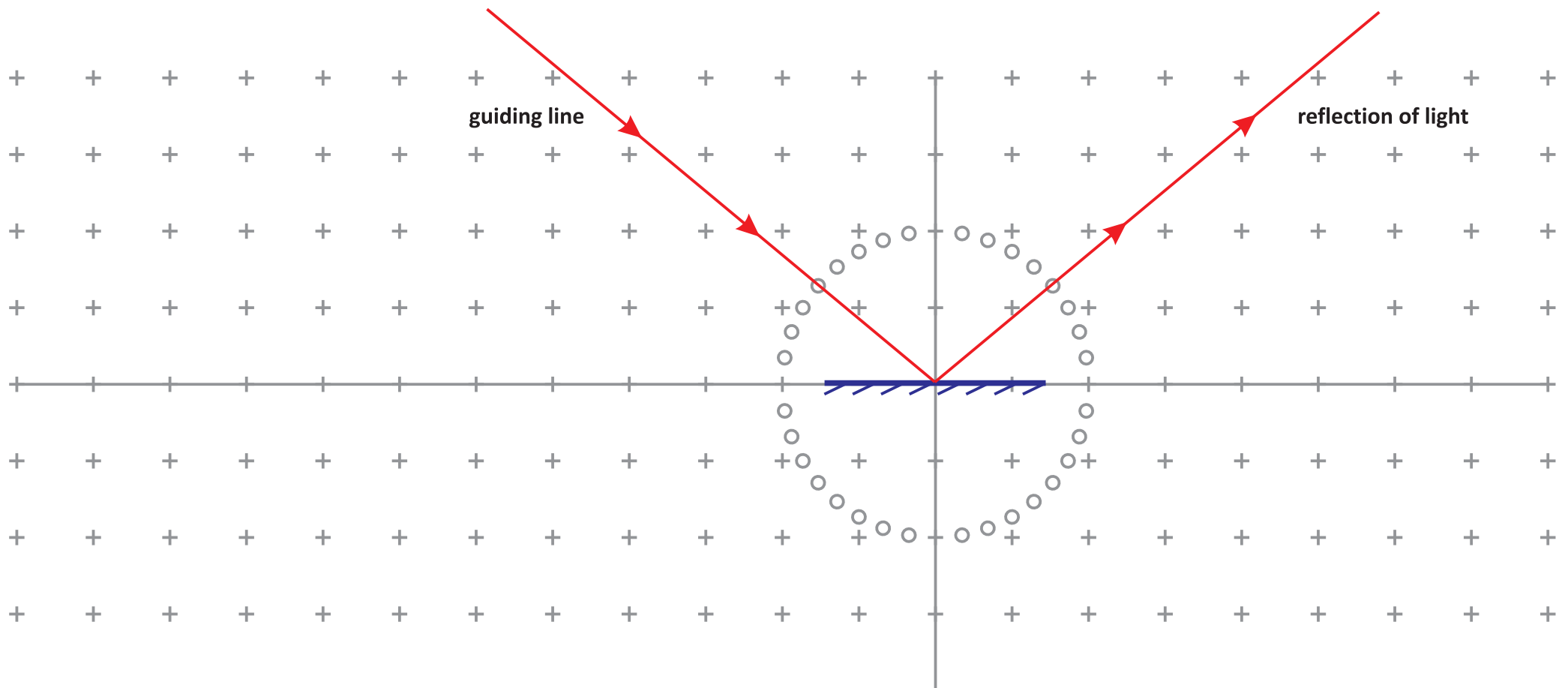

1- Plane Mirror

2- Rotation of mirror and reflection angle


- Place the plane mirror on the determined position in the figure.
- Place the light source and illuminate the plane mirror along the **guiding line**. Keep the incident ray fixed.
- How much is the change in the reflection angle, when you rotate the mirror about the center of protractor (for example by 10 degrees)?

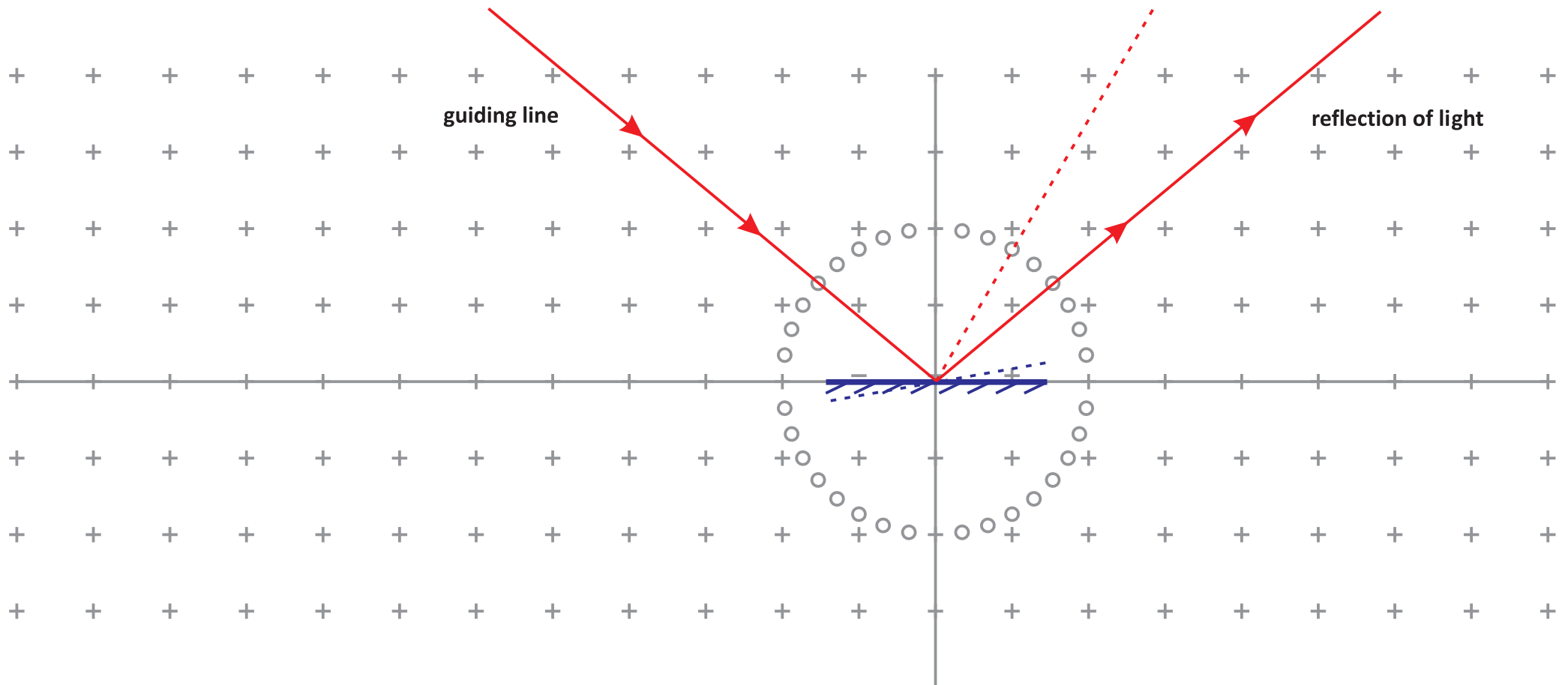

1- Plane Mirror

3- Perpendicular Mirrors


- Place two plane mirrors on the determined position in the figure.
- Illuminate the first mirror along the **guiding line**.
- Reflection ray from the first mirror is the incident ray for the second mirror. What is the situation of the second reflection with respect to the first incident ray?
- Examine different incident angles. Do you know an application for this experiment?

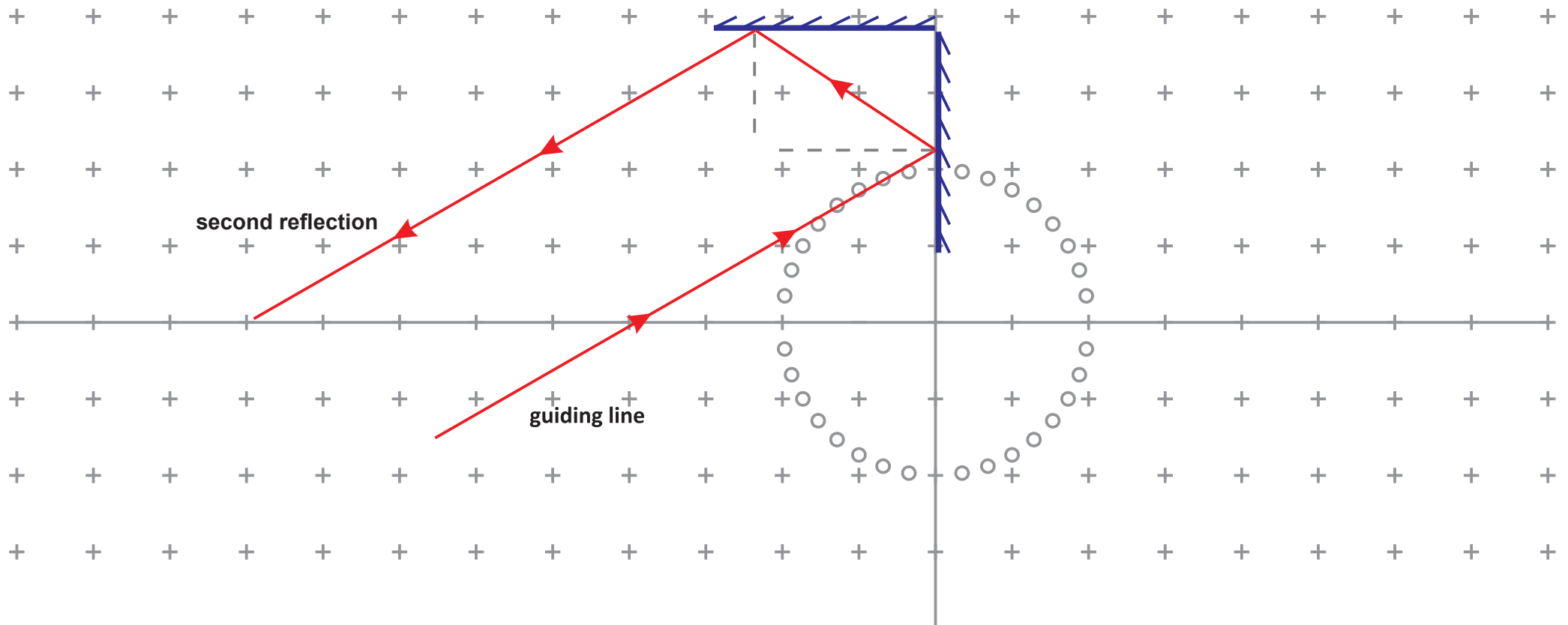

1- Plane Mirror

4- Parallel Mirrors


- Place two plane mirrors on the determined position in the figure.
- Throw the light ray on the first mirror along the guiding line.
- Reflection ray from the first mirror is the incident ray on the second mirror. What is the situation of the second reflection with respect to the first incident ray?
- Repeat the experiment with different incident angles. Do you know an application for this experiment?

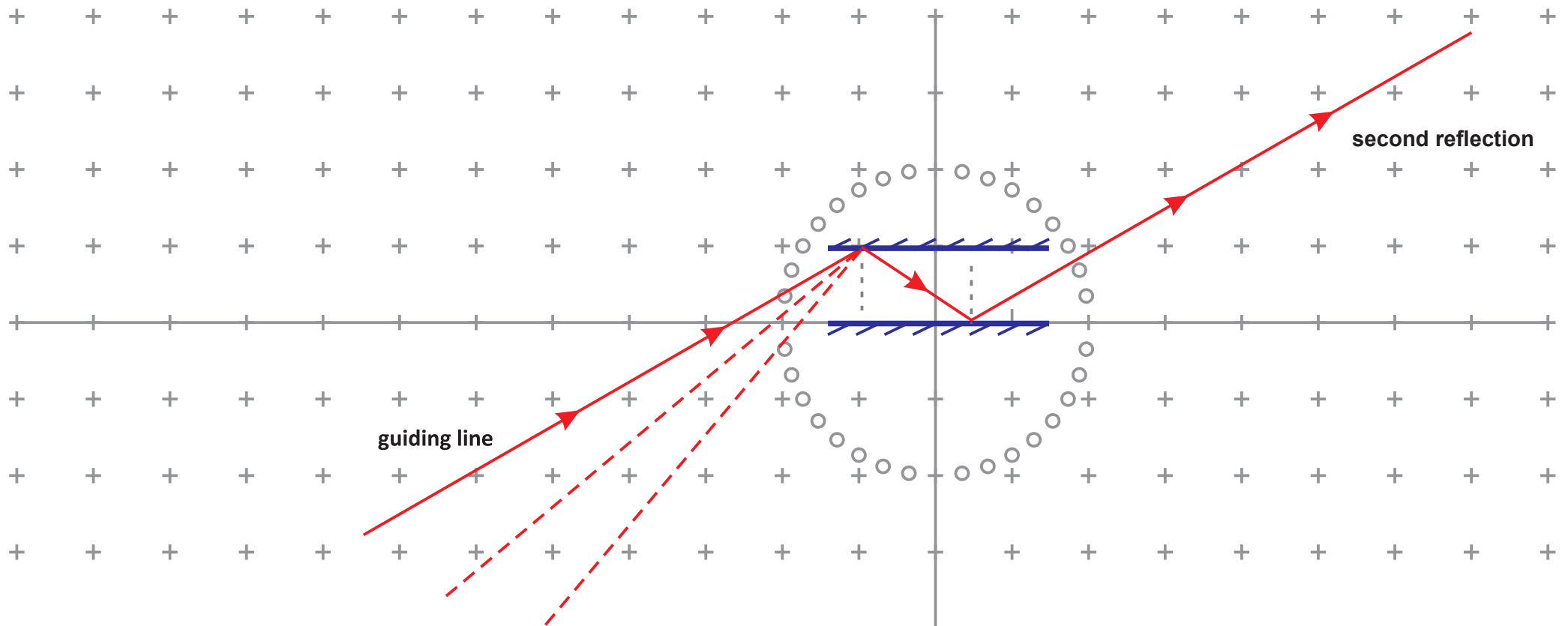

1- Plane Mirror

5- Periscope (Submarine Eye)


- Place two plane mirrors on the determined position in the figure.
- Throw two light rays on the first mirror along the guiding lines.
- Reflection rays from the first mirror are the incident rays on the second mirror. What is the situation of the second reflection with respect to the first incident rays?
- Can you find an application for this experiment?

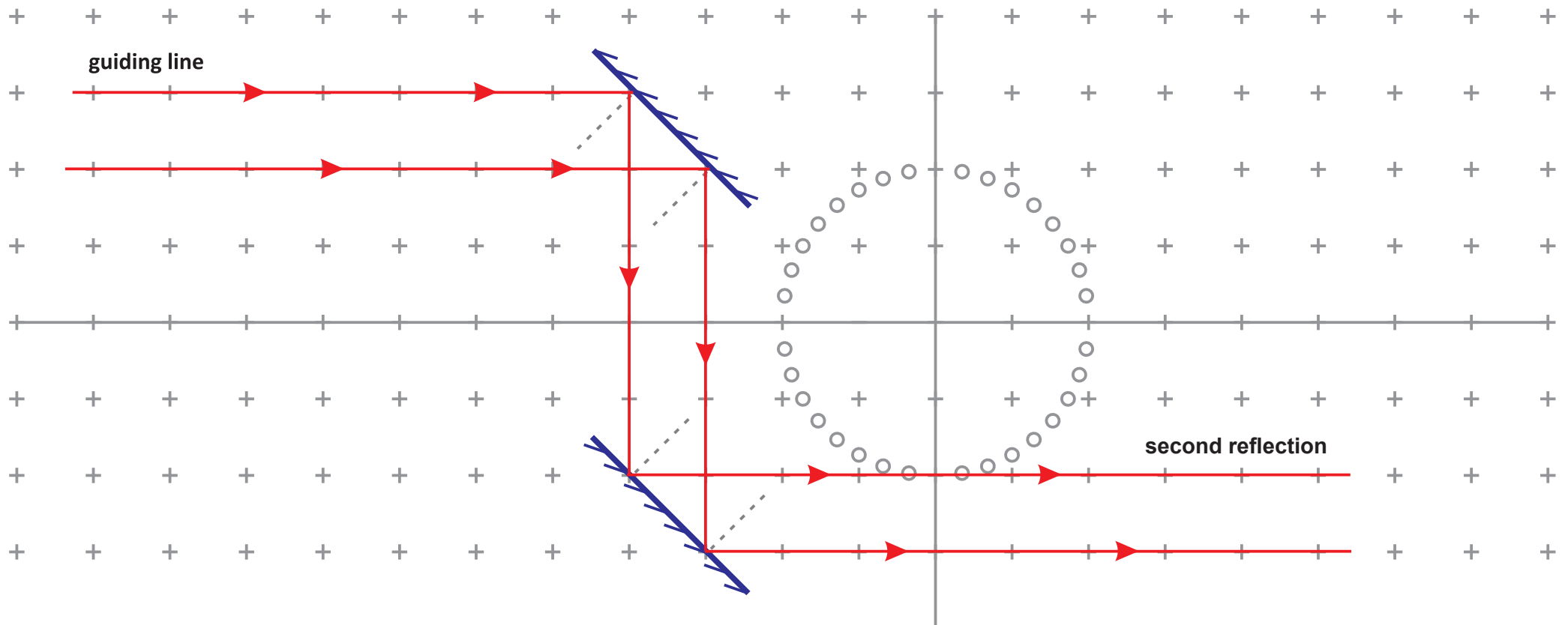

1- Plane Mirror

6- Virtual Image from Real Source


- Place a plane mirror on the determined position in the figure.
- Light rays pass through the point A1 and fall on the mirror. We assume A1 as the real source of rays.
- Reflection rays from the mirror diverge such that it seems they are emanating from the point A2.
- In this case we name A1 and A2, real source and its virtual image, respectively. Why?

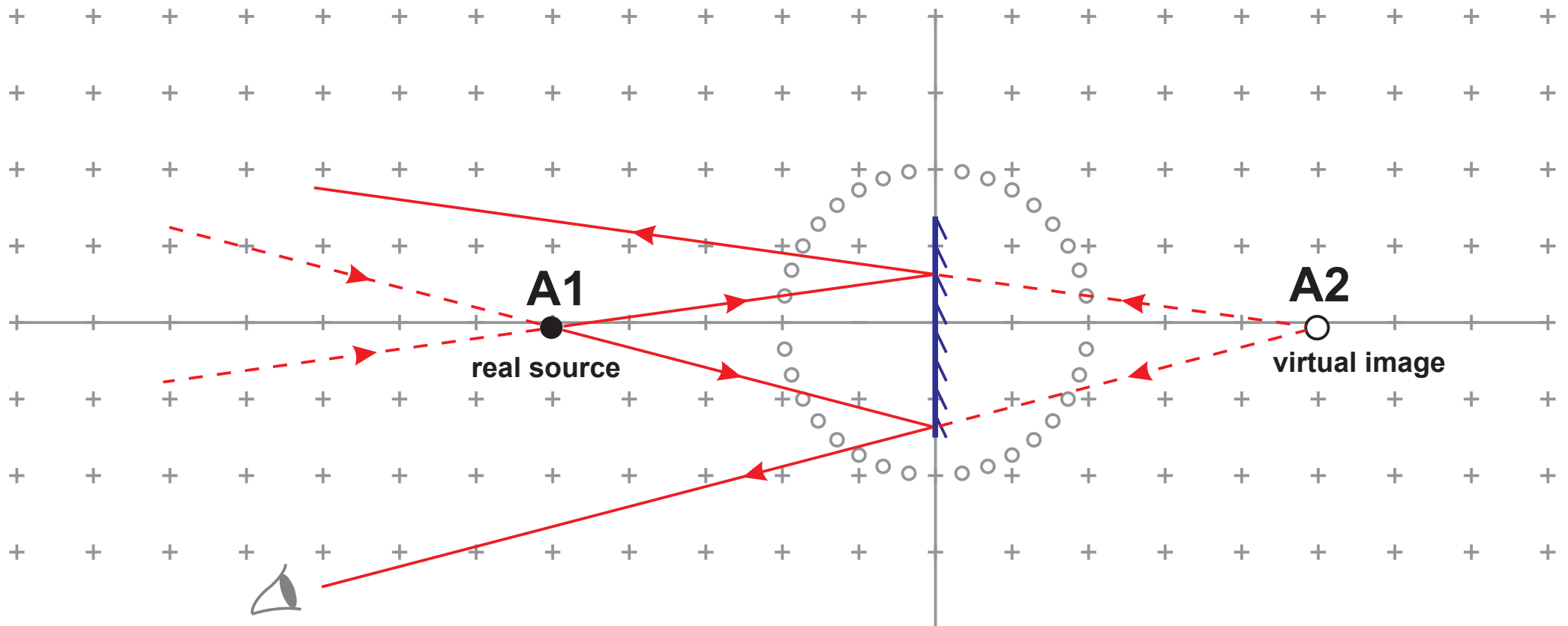

1- Plane Mirror

7- Virtual Image (Hypothetical light source behind the mirror)


- Place the plane mirror on the determined position. Place the mirrorlet (small mirror with a handle) on the point A1.
- Light ray falls on mirrorlet at the point A1. Reflection rays from the mirrorlet illuminate the plane mirror.
- Oscillate the mirrorlet about an axis perpendicular to the board to make incident rays fall on and reflect from the mirror with different angles.
- Reflection rays from the mirror diverge such that it seems they are emanating from the point A2.

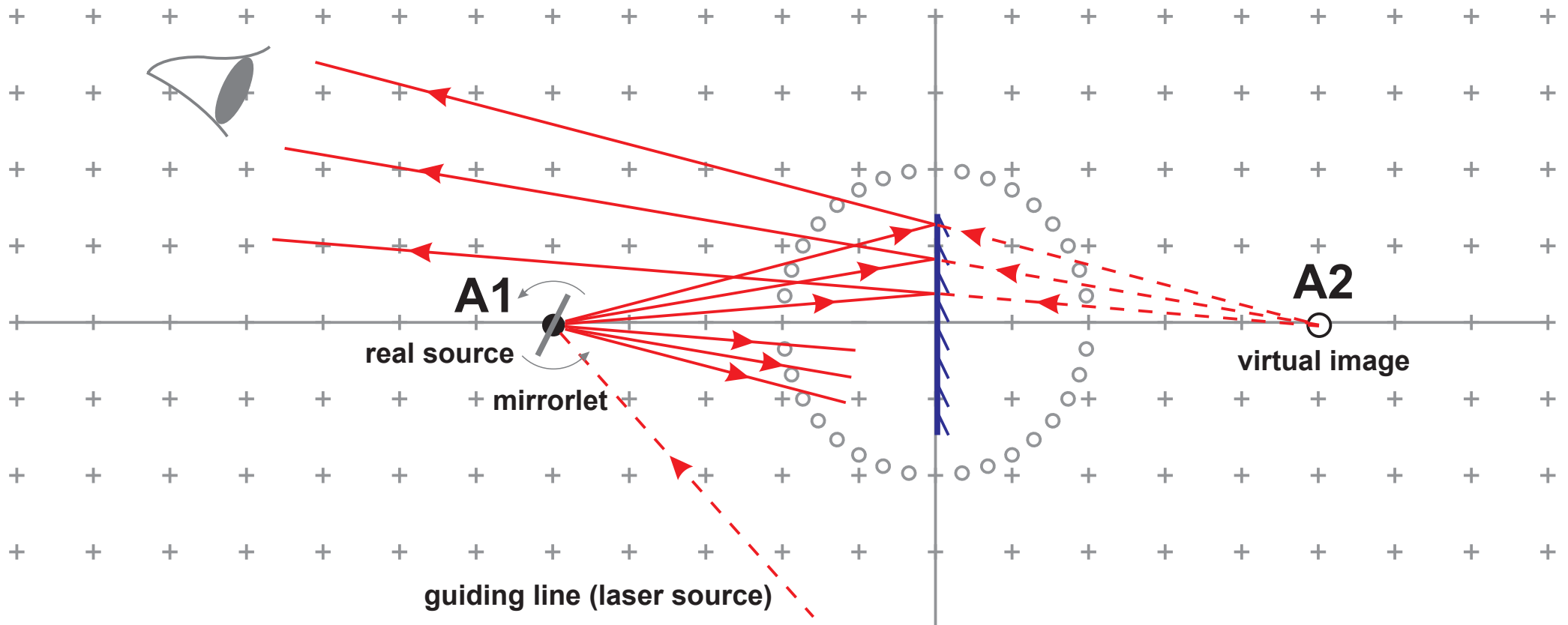

2- Concave Mirror

1- Mirror Vertex


- Place the concave mirror on the determined position in the figure.
- Place a light source on the board such that the incident ray falls on the mirror vertex.
- Make sure to set the mirror position precisely and set the incident ray to fall on the vertex.
- It is seen that the **incident** ray falling on the vertex and its **reflection**, are **symmetric** with respect to the principal axis of concave mirror.

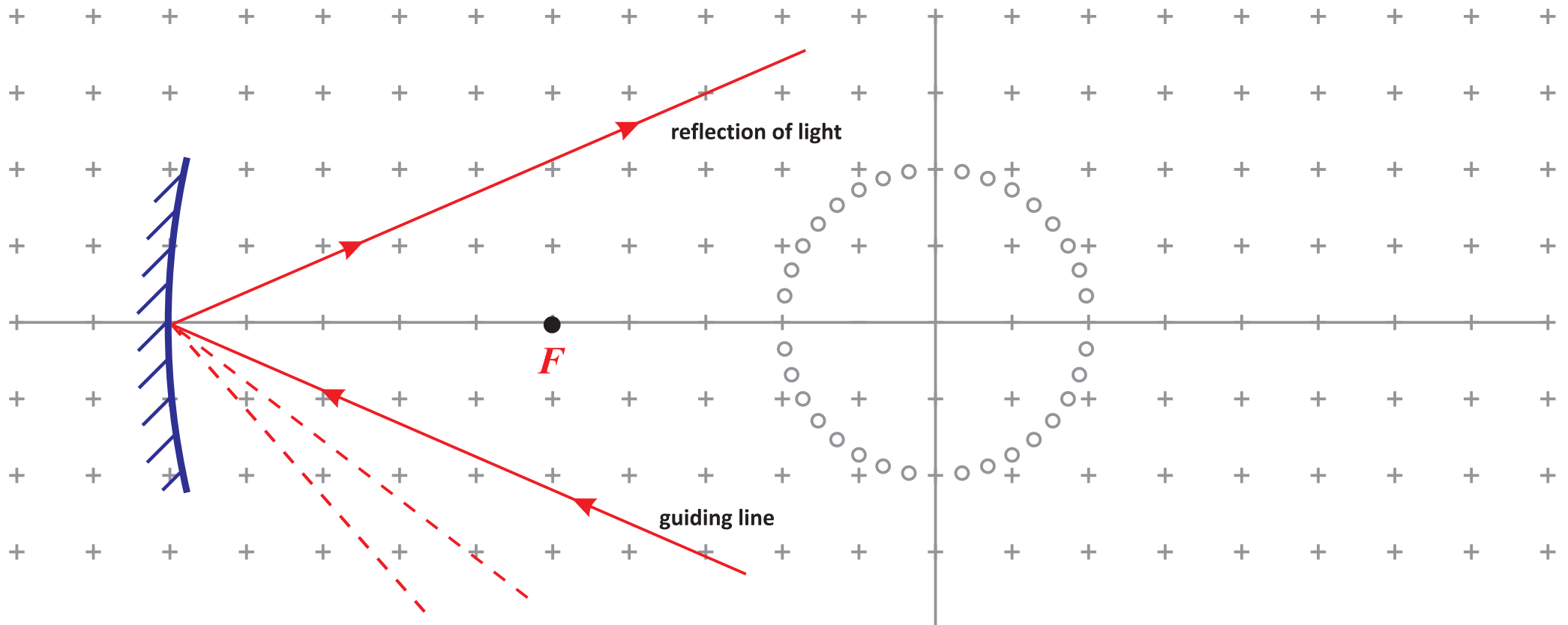

2- Concave Mirror

2- Mirror Center


- Place the concave mirror on the determined position in the figure. Place two or three light sources on the board.
- Illuminate the mirror such that the incident rays pass the mirror center (C) before falling on mirror.
- Make sure to set the mirror position precisely, and that incident rays pass the center.
- It is seen that reflection of the incident rays passing through the center, will coincide with the incident ray.

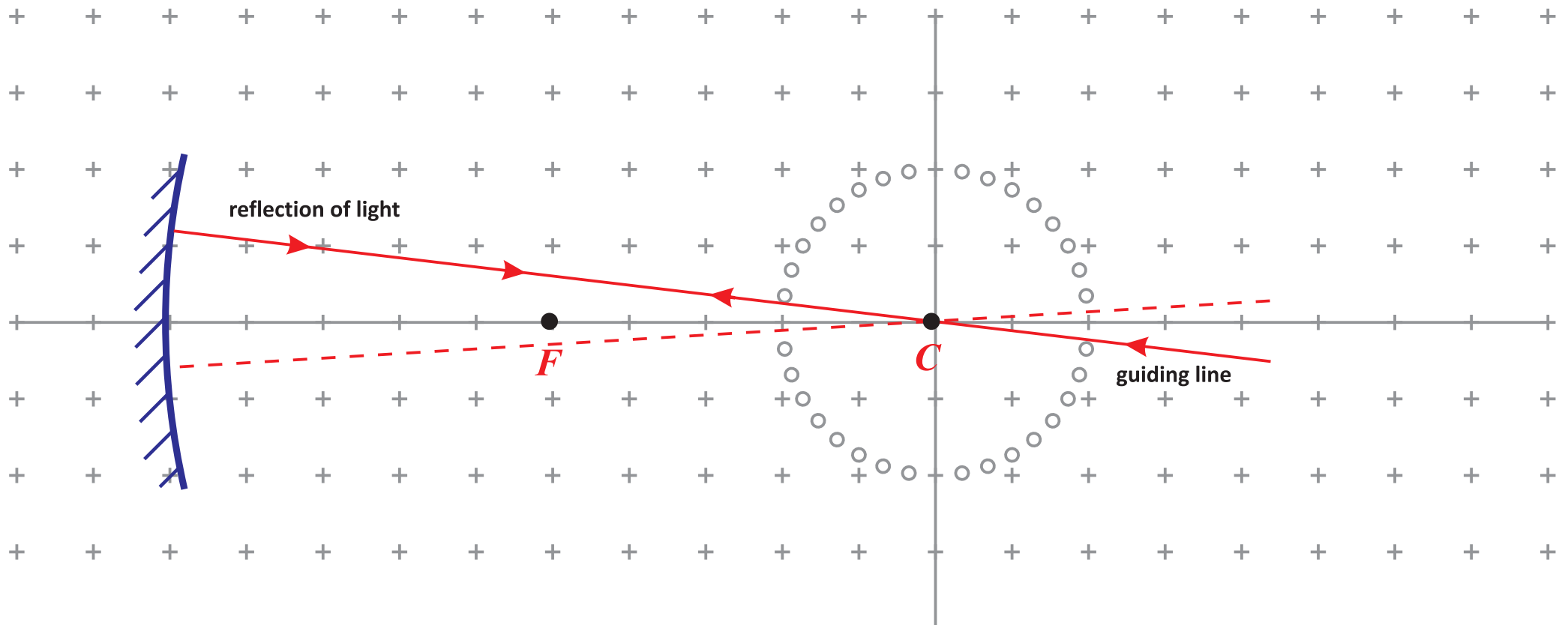

2- Concave Mirror

3- Principal Focal point (Reflection from the incident rays parallel to the principal axis pass through this point)


- Place the concave mirror on the determined position in the figure. Place two or three light sources on the board.
- Illuminate the mirror, by the light rays **parallel to the principal axis** (for example, along the guiding lines).
- Make sure to set the mirror position precisely and set the incident rays parallel to principal axis.
- Reflections of the incident rays parallel to the axis pass through a common point known as the **focal point** of concave mirror (F).

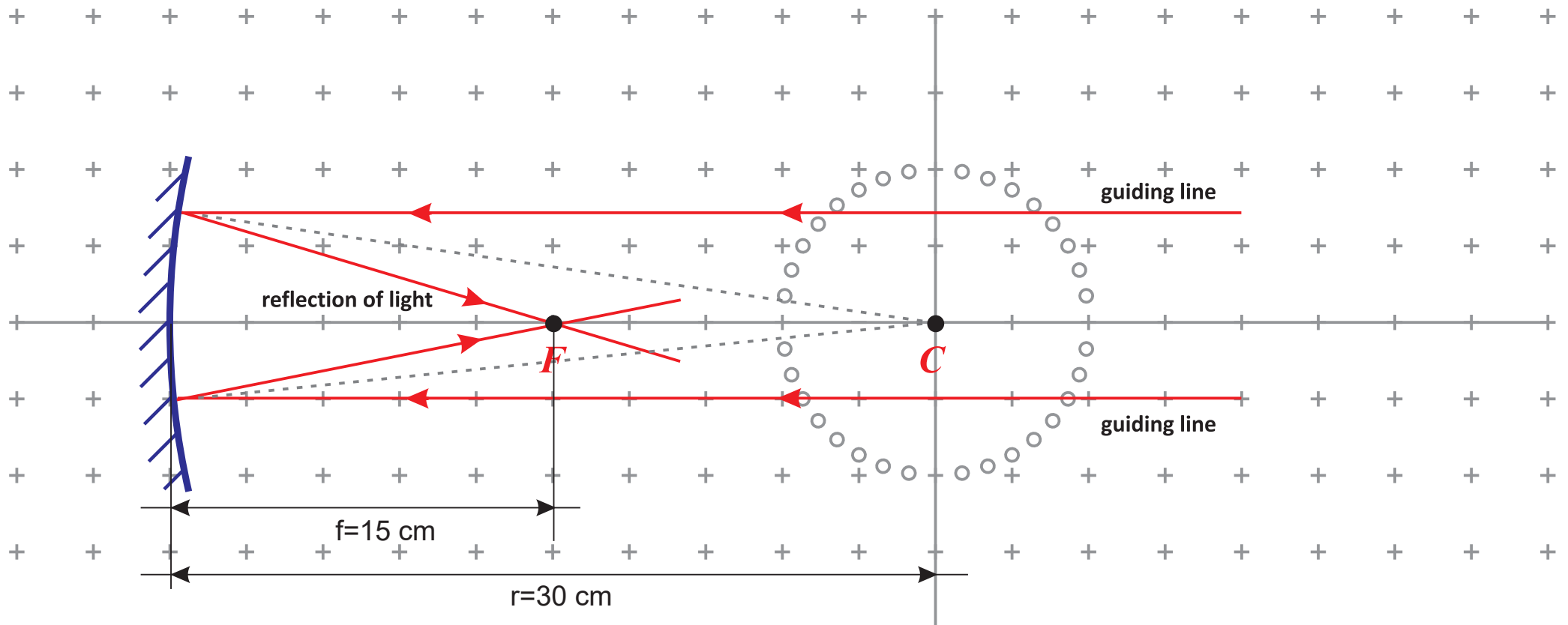

2- Concave Mirror

4- Real Image (Magnification equal to unity)

- Place the concave mirror on the determined position in the figure. Place two or three light sources on board.
- Incident rays pass **A1** and then fall on the mirror. We assume **A1** to be the **real source** of rays.
- Reflections converge at the point **A2** and **A2** becomes a new light source.
- In this case we call **A1** a **real source** and **A2** its **real image**. Why?

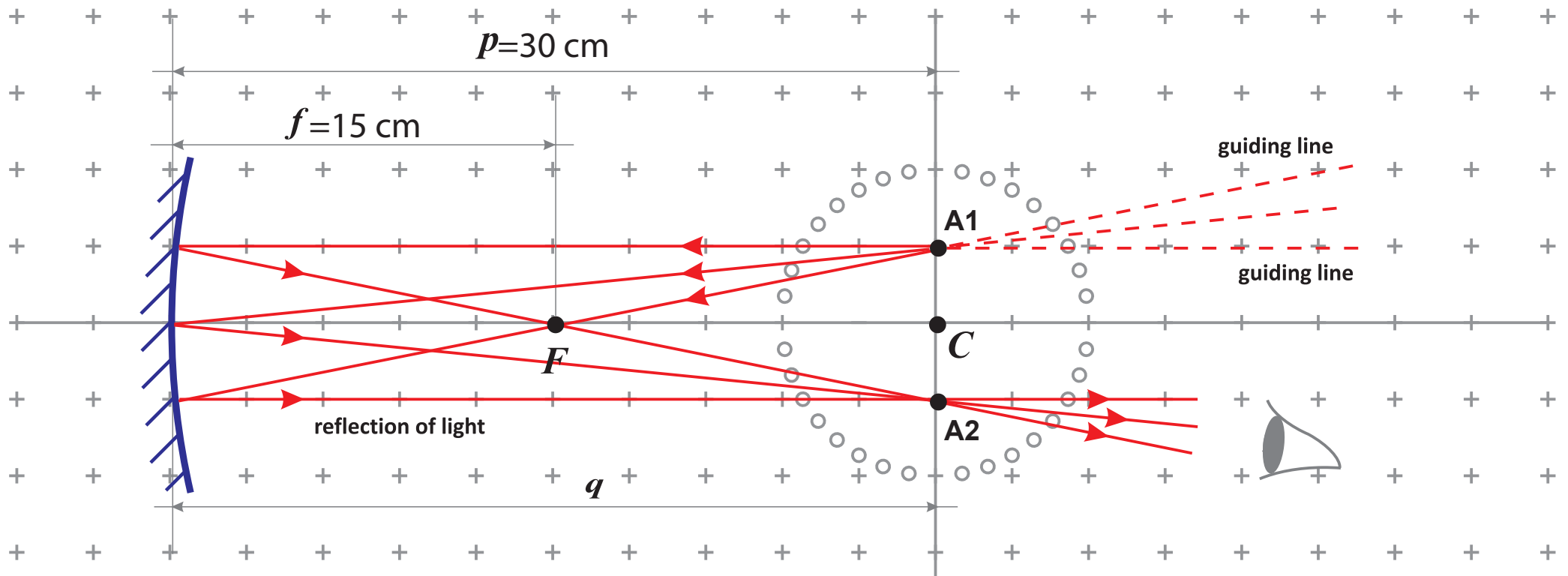


$$\frac{1}{p} + \frac{1}{q} = \frac{1}{f} \Rightarrow \frac{1}{30 \text{ cm}} + \frac{1}{q} = \frac{1}{15 \text{ cm}} \Rightarrow q = 30 \text{ cm}$$

2- Concave Mirror

5- Real Image (Object: Infinite number of rays, Image: Infinite number of rays)

- Place a concave mirror on the determined position in the figure. Place a mirrorlet on the point A1.
- Source light ray falls on mirrorlet at the point A1. Reflection rays from the mirrorlet illuminate the concave mirror.
- Oscillate the mirrorlet about an axis perpendicular to the board to make incident rays fall on, and reflect from the mirror with different angles.
- A2 is the real image of A1. From each point of source and its image (real or virtual), infinite number of rays are emitted.

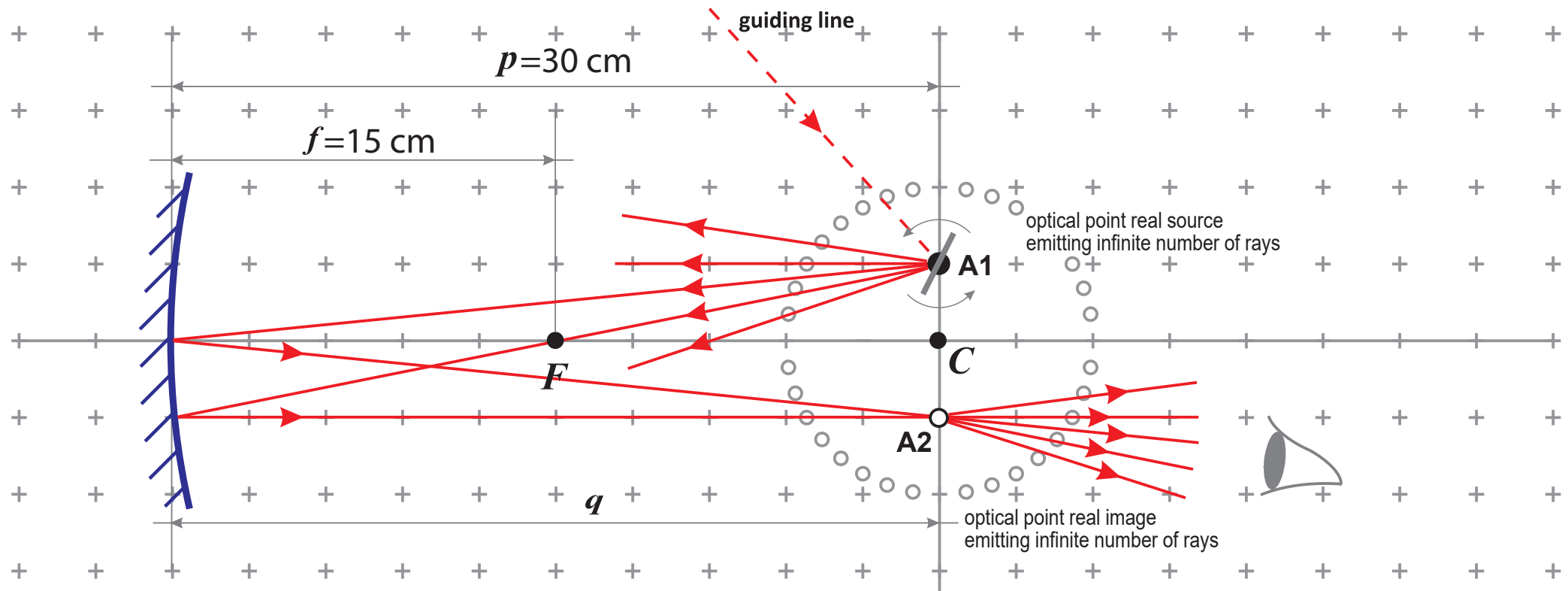


$$\frac{1}{p} + \frac{1}{q} = \frac{1}{f} \Rightarrow \frac{1}{30\text{ cm}} + \frac{1}{q} = \frac{1}{15\text{ cm}} \Rightarrow q = 30\text{ cm}$$

2- Concave Mirror

6- Real Image (Object: Infinite number of sources, Each source: Infinite number of rays)

- Place a plane mirror on the determined position in the figure. Place the mirrorlet on the point A1.
- By oscillating the mirrorlet emit light rays from A1 to the concave mirror and get their reflection.
- A2 is the real image of A1. Repeat this experiment for other points of the optical object A1B1C1D1.
- Each optical object and its image are consisted of infinite number of point sources and from each source infinite number of rays is emitted.

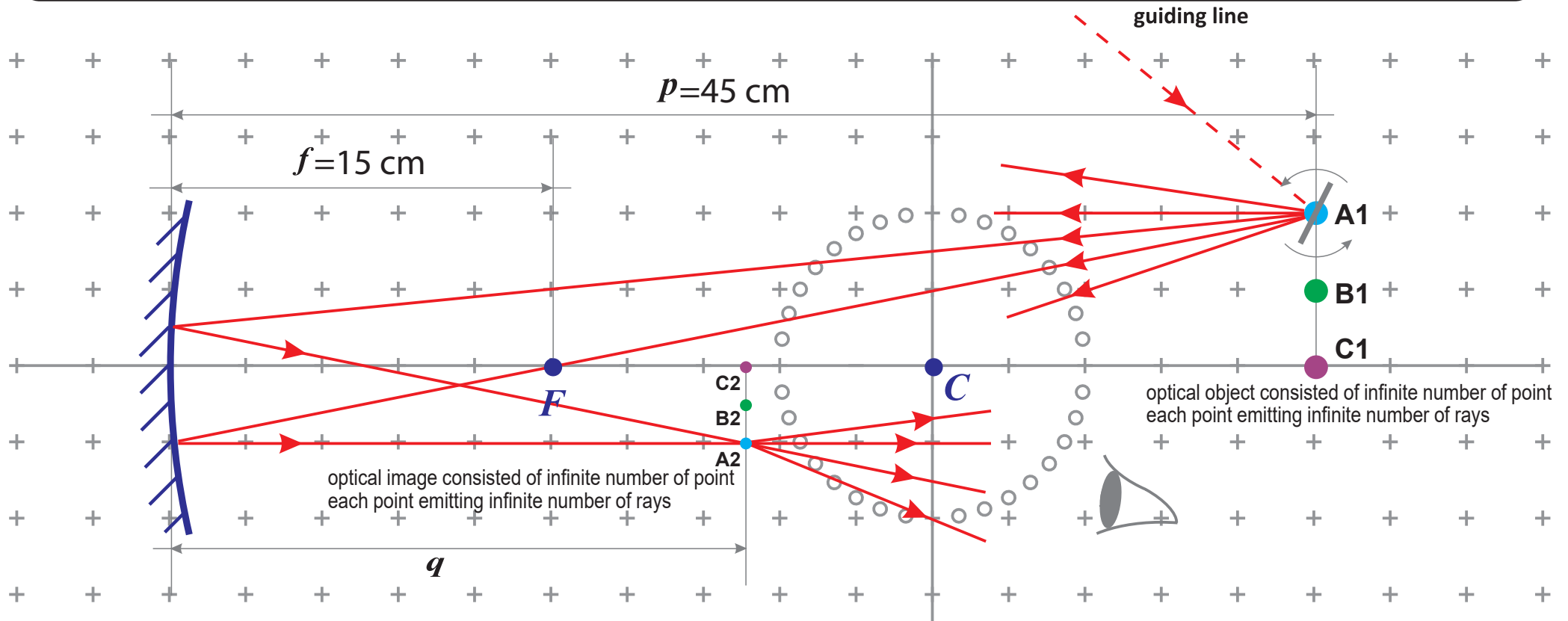


$$\frac{1}{p} + \frac{1}{q} = \frac{1}{f} \Rightarrow \frac{1}{45 \text{ cm}} + \frac{1}{q} = \frac{1}{15 \text{ cm}} \Rightarrow q = 22.5 \text{ cm}$$

3- Convex Mirror

1- Mirror Vertex


- Place the convex mirror on the determined position in the figure.
- Place a light source on the board such that the incident ray falls on the mirror vertex.
- Make sure to set the mirror position precisely and set the incident ray to fall on the vertex.
- It is seen that the **incident** ray falling on the vertex and its **reflection**, are **symmetric** with respect to the principal axis of convex mirror.

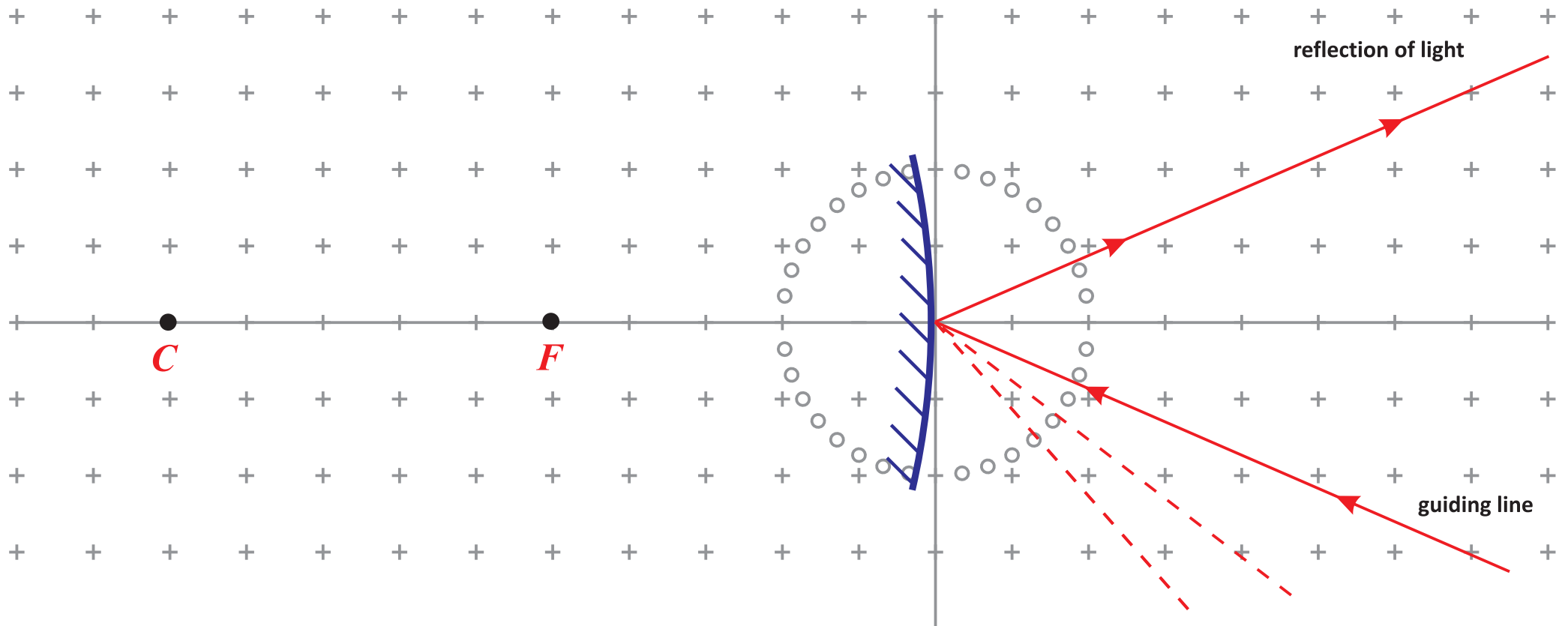

3- Convex Mirror

2- Mirror Center


- Place the convex mirror on the determined position in the figure. Place a light sources on the board.
- With the aid of guiding lines illuminate the mirror such that the incident light rays are directed toward the center (C).
- Make sure to set the mirror position precisely, and that incident ray is directed toward the center.
- It is seen that in the convex mirror, incident ray directing toward the center, reflect on itself.

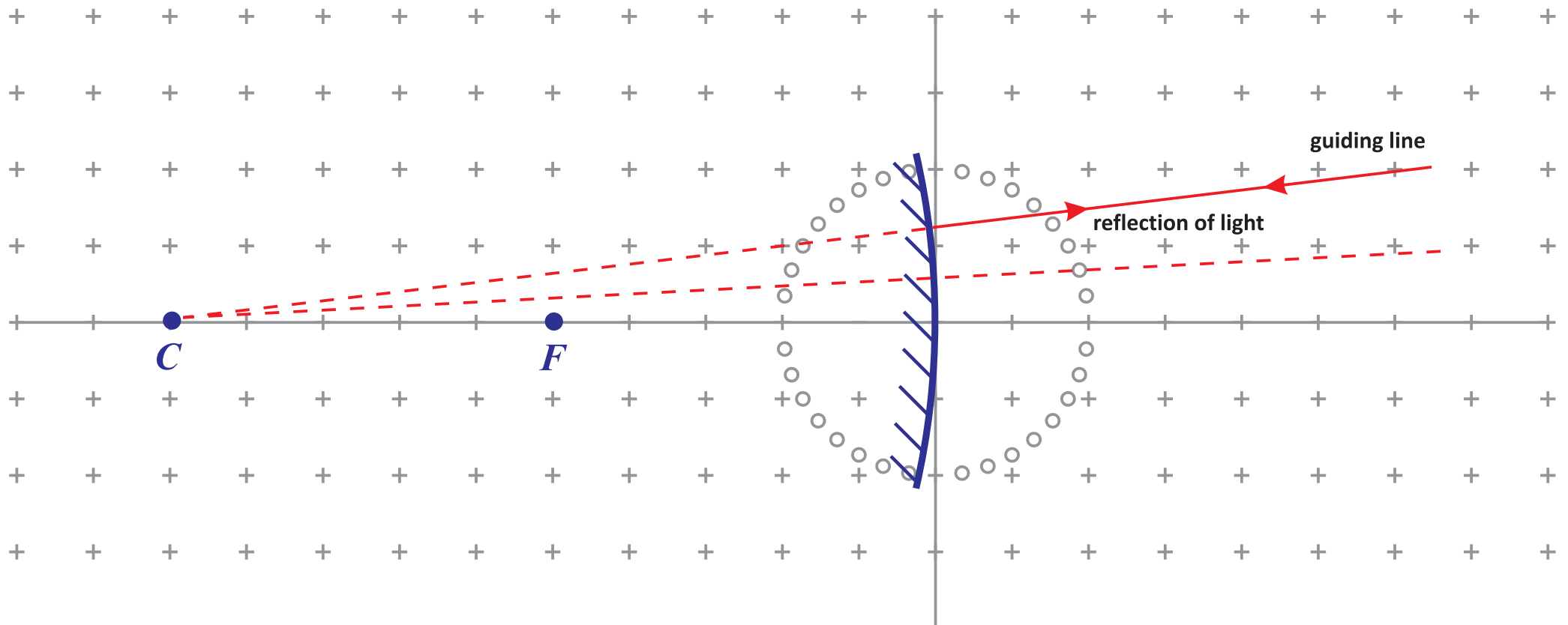

3- Convex Mirror

3- Principal Focal Point (Intersection of the direction of reflections corresponding to incident rays parallel to the principal axis)


- Place the concave mirror on the determined position in the figure. Place two or three light sources on the board.
- Illuminate the mirror, by the light rays **parallel to the principal axis** (for example, along the guiding lines).
- If the direction of reflection of an incident ray parallel to the axis is extended, it will pass through the **focal point** of concave mirror (**F**).

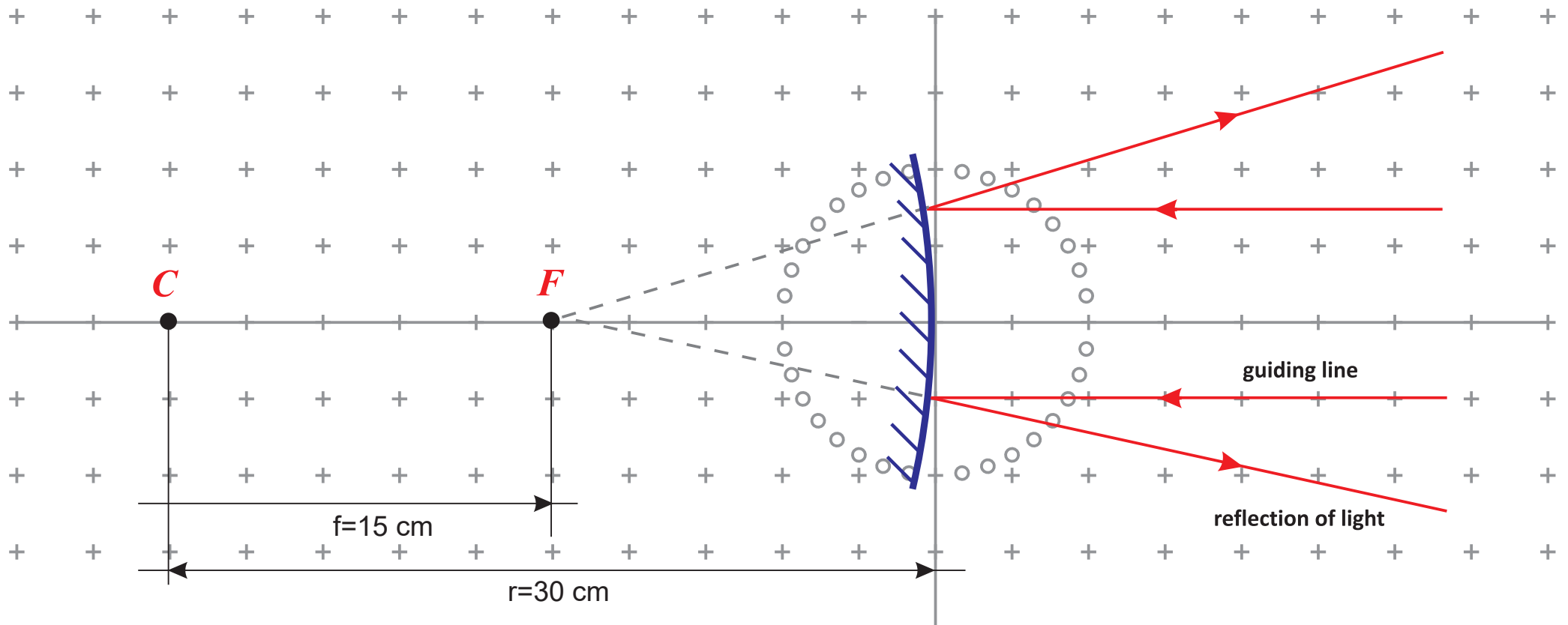

3- Convex Mirror

4- Virtual Image (Magnification less than unity)

- Place the convex mirror on the determined position in the figure. Place two or three light sources on the board.
- Incident rays pass **A1** and fall on the mirror. We assume **A1** to be the **real source** of rays.
- Reflections diverge from the mirror such that it **seems** they are sourced from **A2**.
- In this case **A1** is named a real source and **A2** is called its **virtual image**. Why?

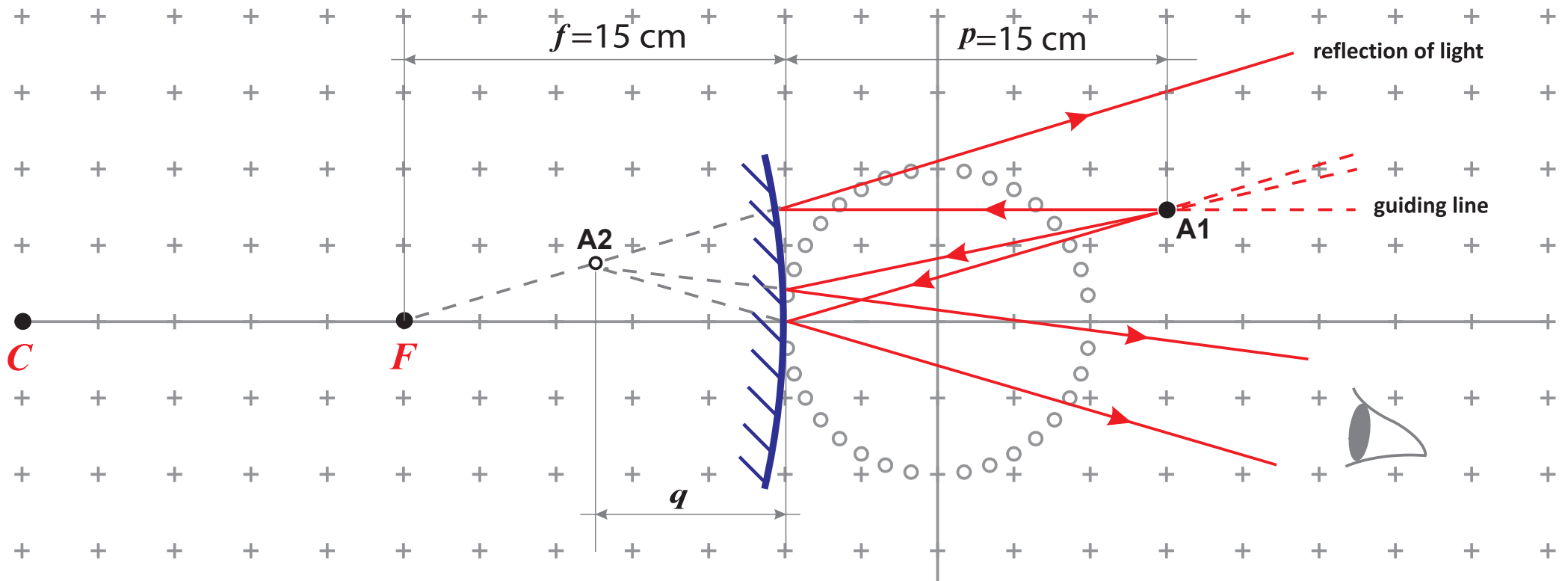


$$\frac{1}{p} + \frac{1}{q} = \frac{1}{f} \Rightarrow \frac{1}{15 \text{ cm}} + \frac{1}{q} = \frac{1}{-15 \text{ cm}} \Rightarrow q = -7.5 \text{ cm}$$

4- Refraction of Light

1- Transmission of light between two transparent media


- Place the transparent half-cylinder on the determined position in the figure.
- Place a light source on the board.
- Throw the light ray on the curved surface of half-cylinder along the radius and toward the center of protractor.
- By moving the light source and changing the **incidence angle**, observe the **refraction** of light when passing from the high density medium to the low density medium (air).

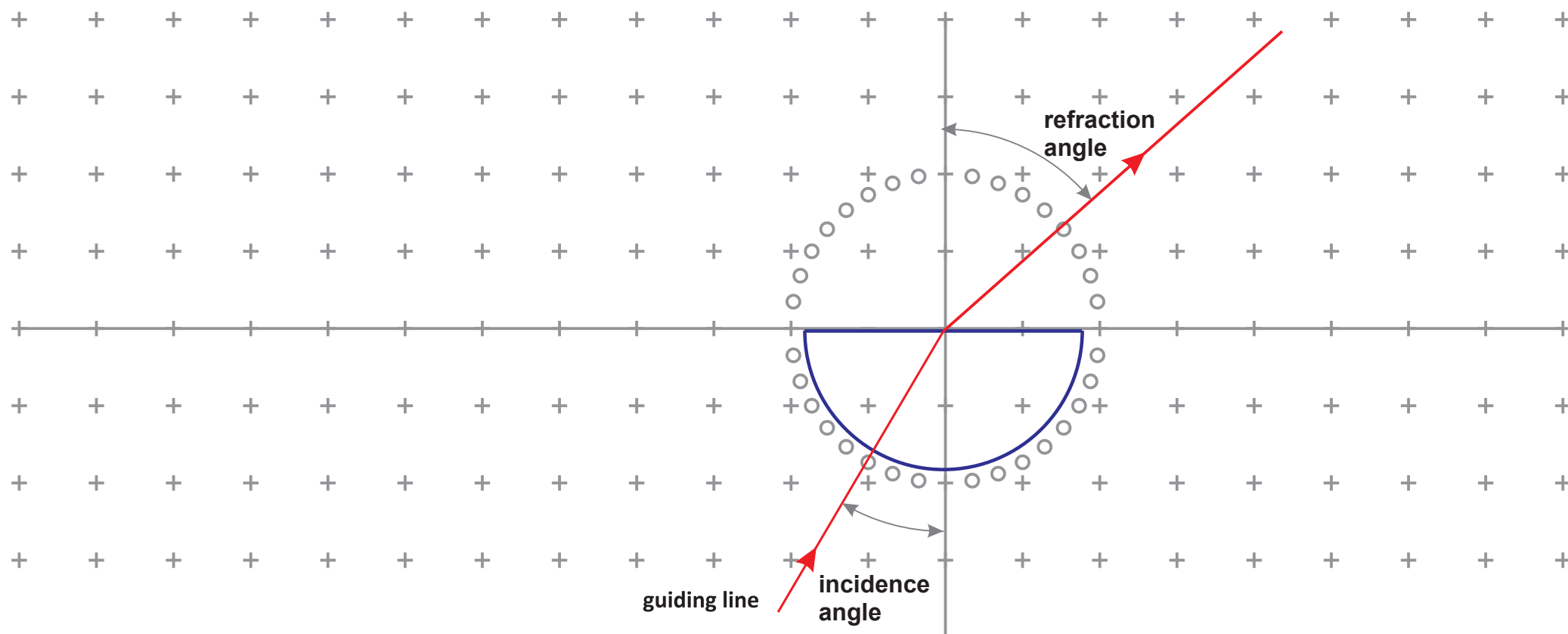

4- Refraction of Light

2- Total reflection, Partial reflection and critical angle


- Place the transparent half-cylinder on the determined position in the figure. Place a light source on the board.
- Throw the light ray on the curved surface of half-cylinder along the radius and toward the the center of protractor.
- By changing the **incidence angle**, for a special incidence angle called the **critical angle**, reflection ray becomes tangent to the flat surface of half-cylinder and for larger incidence angles, total reflection will begin.

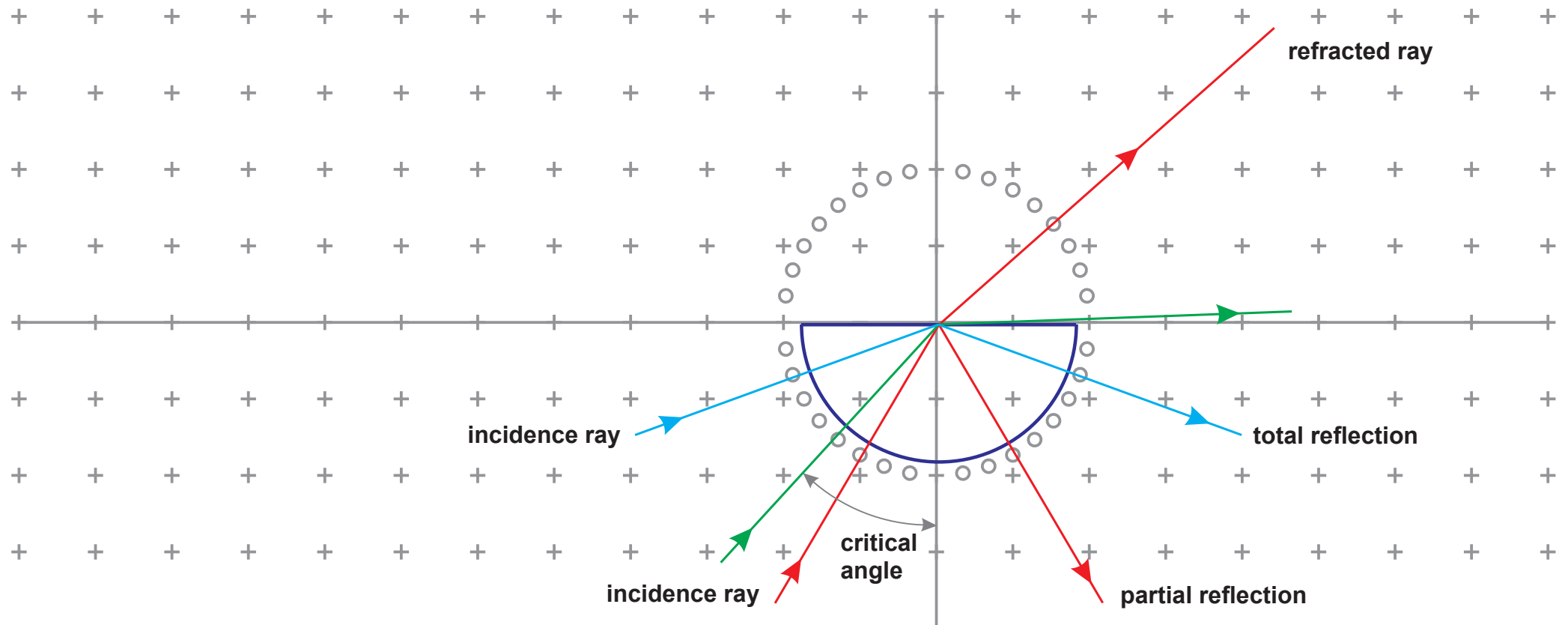

5- Refraction of Light

1- Transmission of light through a parallel-plane slab


- Place the transparent parallel-plane slab on the determined position in the figure.
- Place two light sources on the left side of the board.
- With the aid of guiding lines, illuminate the surface of plane-parallel slab.
- Change the incidence angle of light by rotating the slab in its place and observe the amount of shift in the direction of light rays.

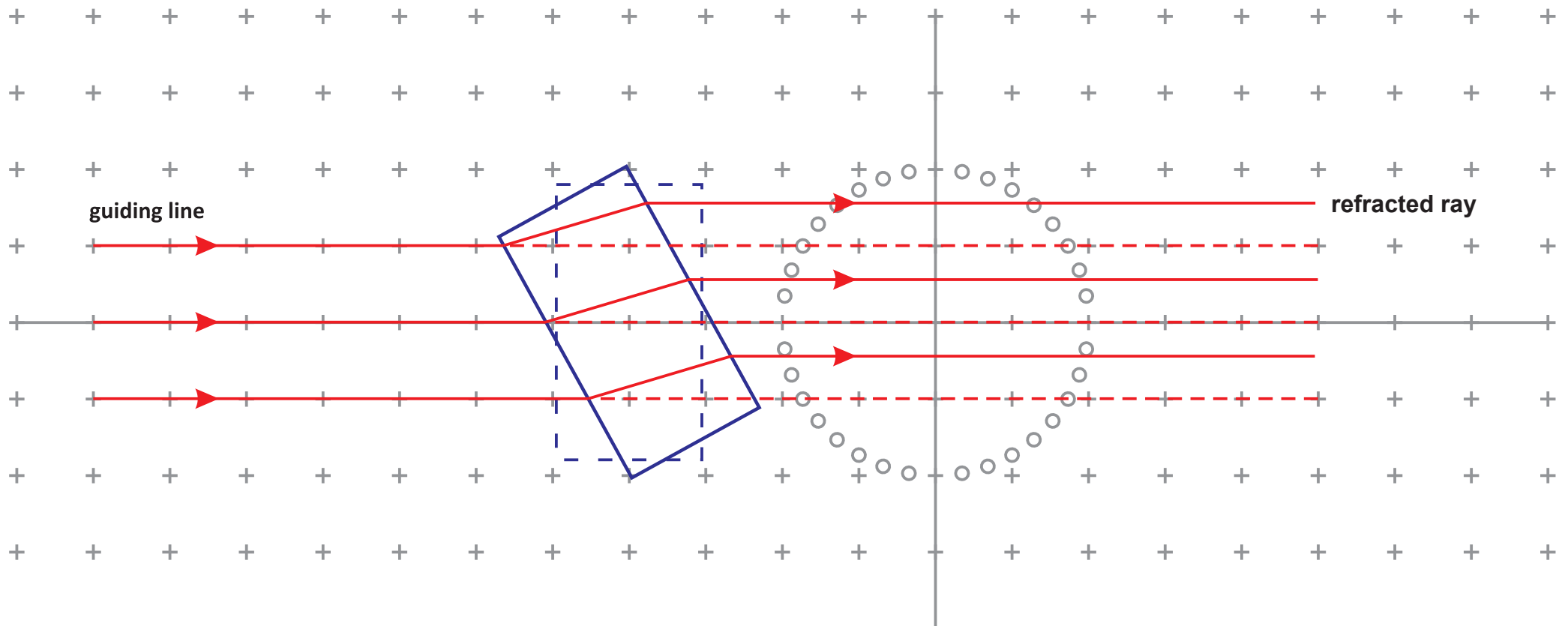

6- Prism

1- Deflection Angle (Refraction of light by prism)


- Place the transparent prism on the determined position in the figure.
- By a light source and with the aid of guiding lines illuminate the surface of prism.
- Observe the refraction of light by prism.
- Angle between the outgoing ray and the direction of ingoing ray is called the light **deflection angle**.

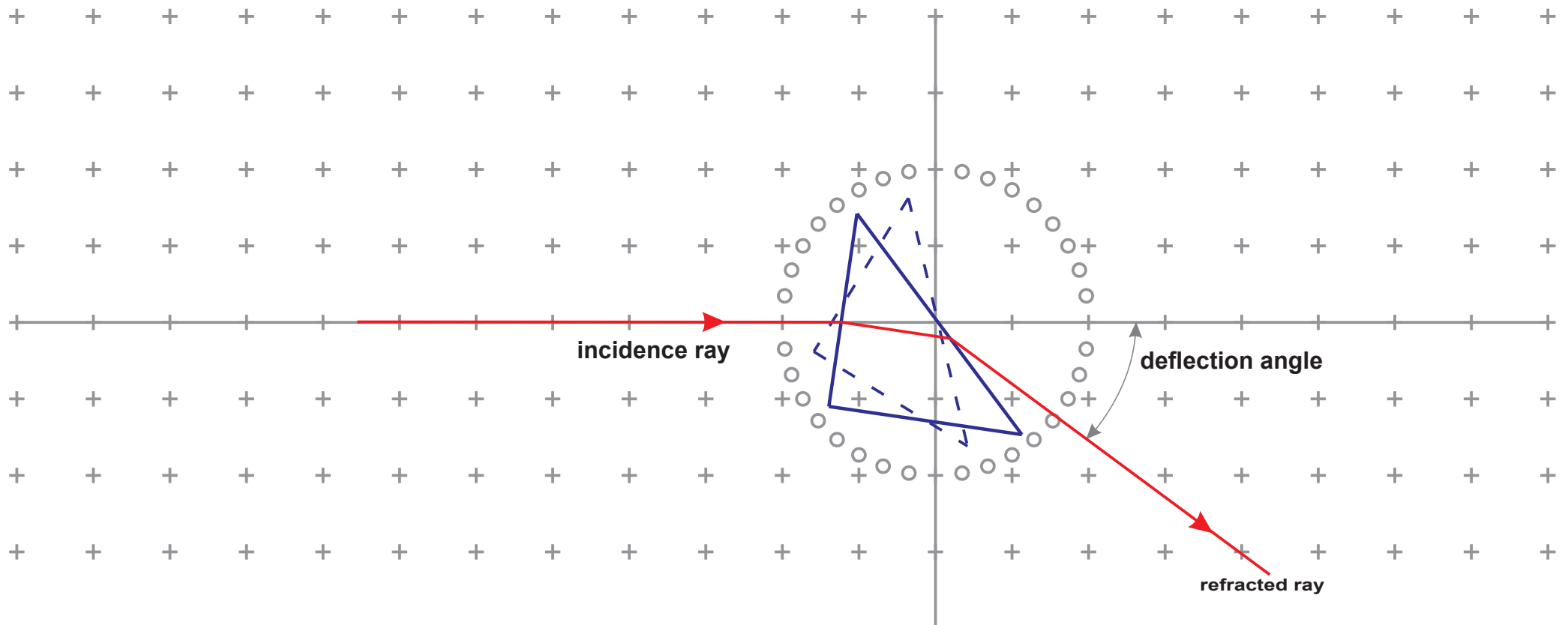

6- Prism

2- A total Reflection (90 degrees light deflection by the prism)


- Place the transparent prism on the determined position in the figure.
- By a light source and with the aid of guiding lines illuminate the prism and observe the result.
- In this case, a **total reflection** occurs instead of refraction, and light is deflected by 90 degrees.
- This phenomenon shows that the **critical angle** for the medium is less than the incidence angle of 45 degrees.

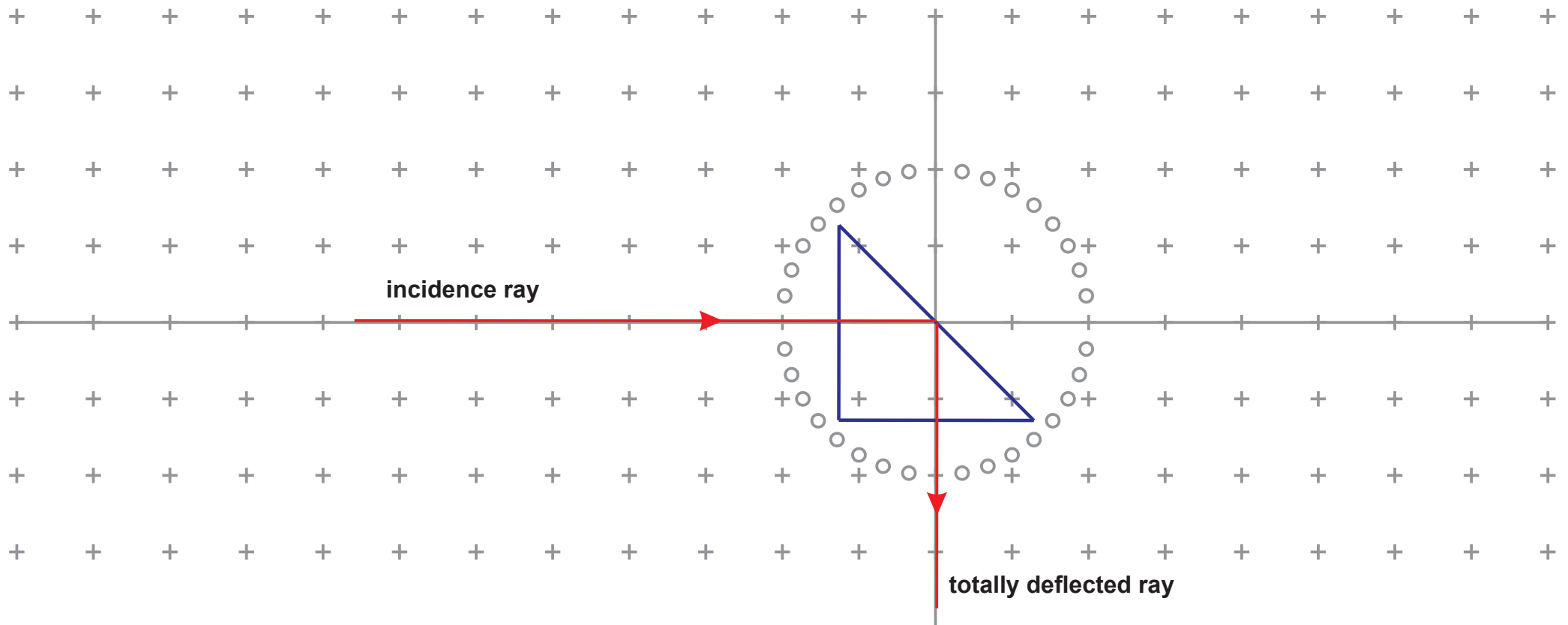

6- Prism

3- Two Total Reflections (180 degrees light deflection by the prism)


- Place the transparent prism on the determined position in the figure.
- By a light source and with the aid of guiding lines illuminate the prism and observe the result.
- In this case, **two total reflections** occur instead of refraction, and light is deflected by 180 degrees.
- This phenomenon shows that the **critical angle** for the medium is less than the incidence angle of 45 degrees.

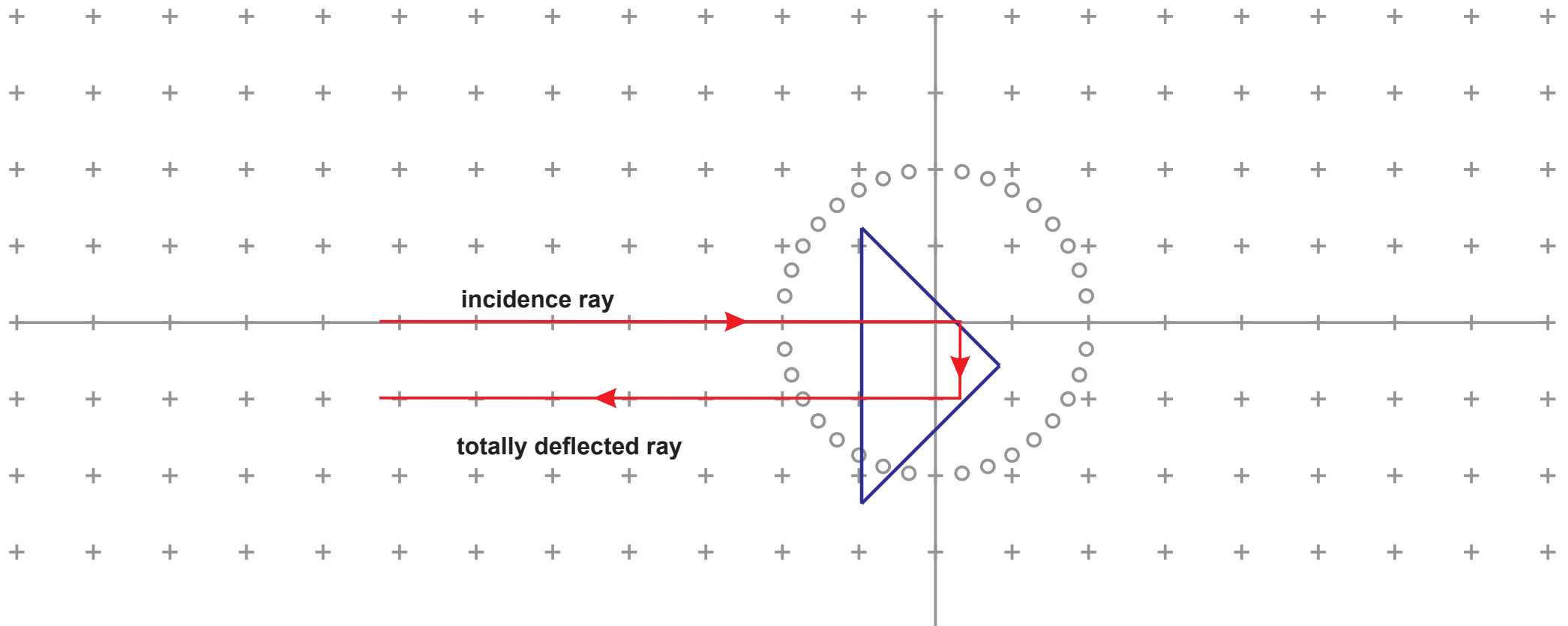

7- Convex Lens

1- Optical Center


- Place a convex lens on the determined position in the figure. Place two or three light sources on the board.
- With the aid of guiding lines set the incident rays such that they fall on the **optical center** of lens.
- Make sure that the lens position is precise, and that the light rays pass through the optical center.
- Each incident ray falling on the optical center, will pass through the lens **without refraction** and deviation in its path.

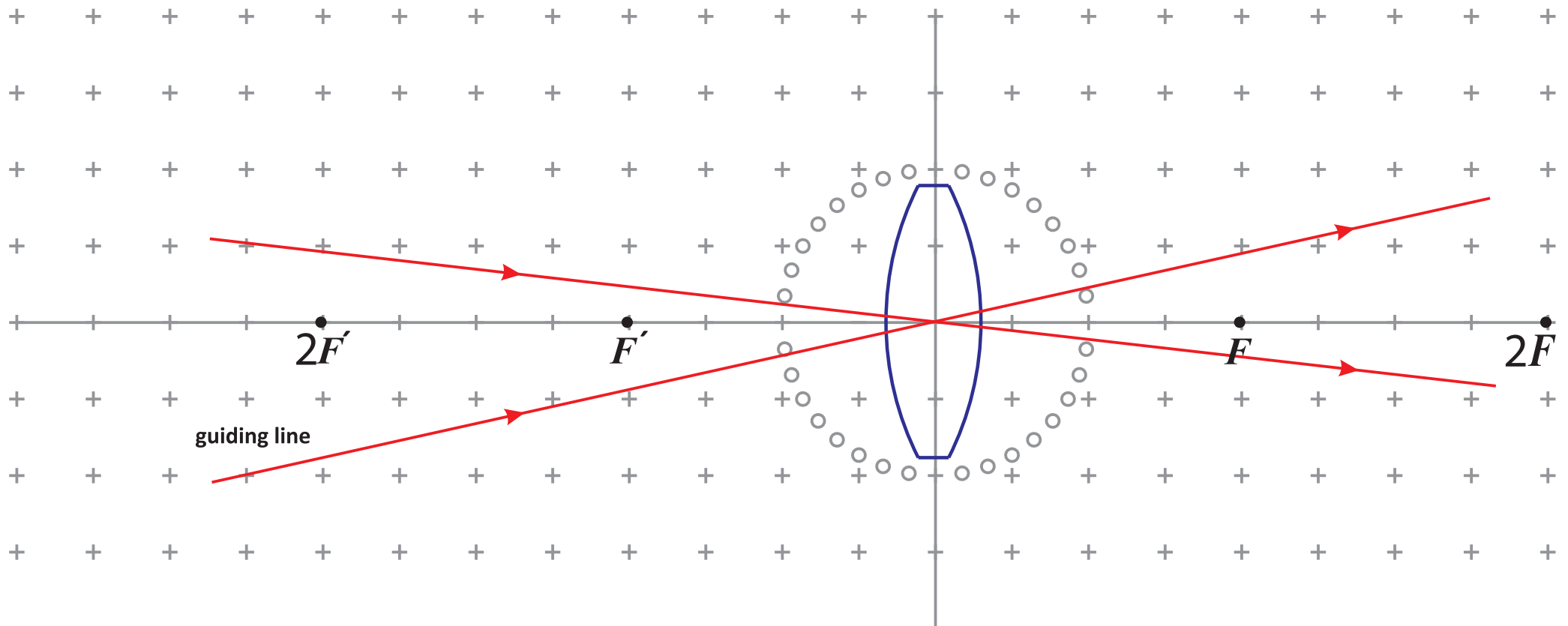

7- Convex Lens

2- Principal Focal Point (Collimation point of the light rays parallel to the principal axis after refraction)


- Place the convex lens on the determined position in the figure.
- by two or three light sources, illuminate the lens with light rays parallel to the principal axis.
- Make sure that the lens position is precise, and that the light rays are parallel with the principal axis.
- Incident rays parallel to the principal axis, after refraction pass through a common point called the convex mirror **focal point (F)**.

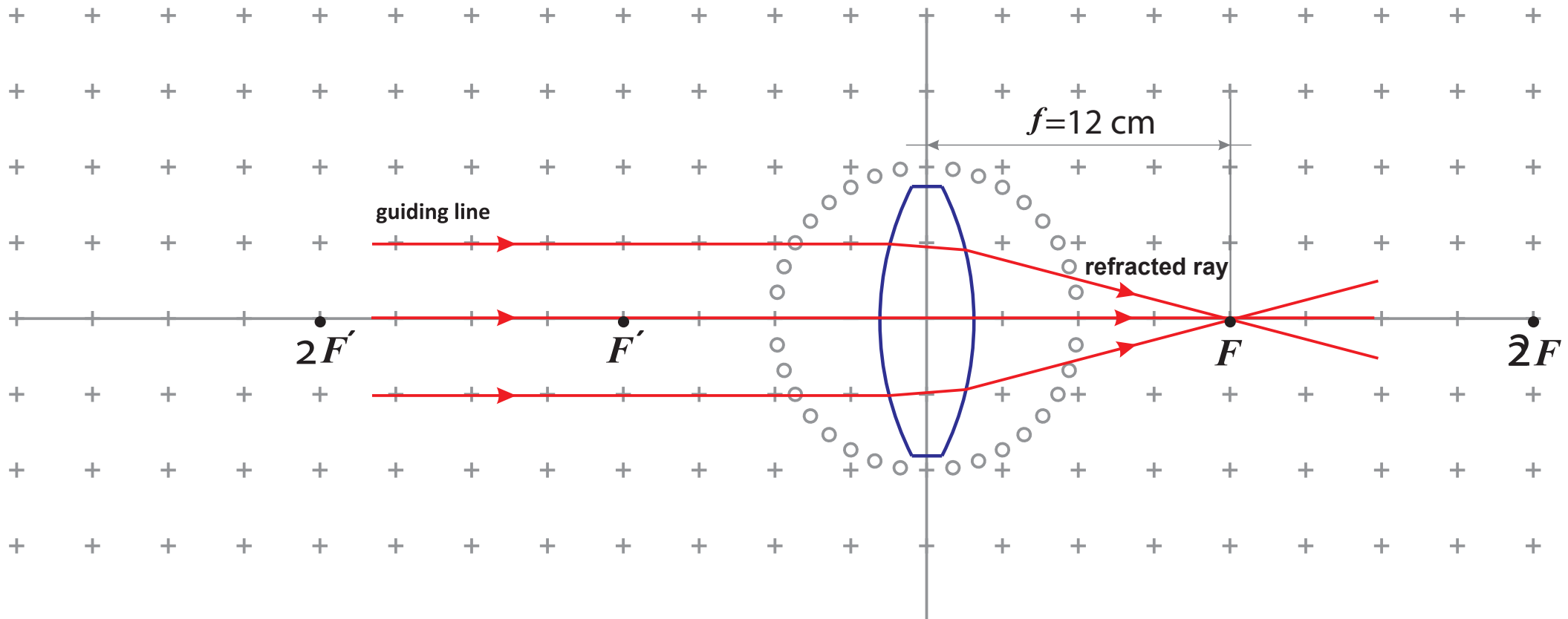

7- Convex Lens

3- Principal Focal Point (Effect of convexity on the position of focal point)


- First place a **plano-convex** lens and then **two plano-convex** lenses on the determined position in the figure. Then by two or three light sources, illuminate the lens with light rays parallel to the principal axis. (Two different experiments)
- Make sure that the lens position is precise, and that the light rays parallel with the principal axis.
- Incident rays parallel to the principal axis, pass through a common point called the convex mirror focal point (F), after refraction.

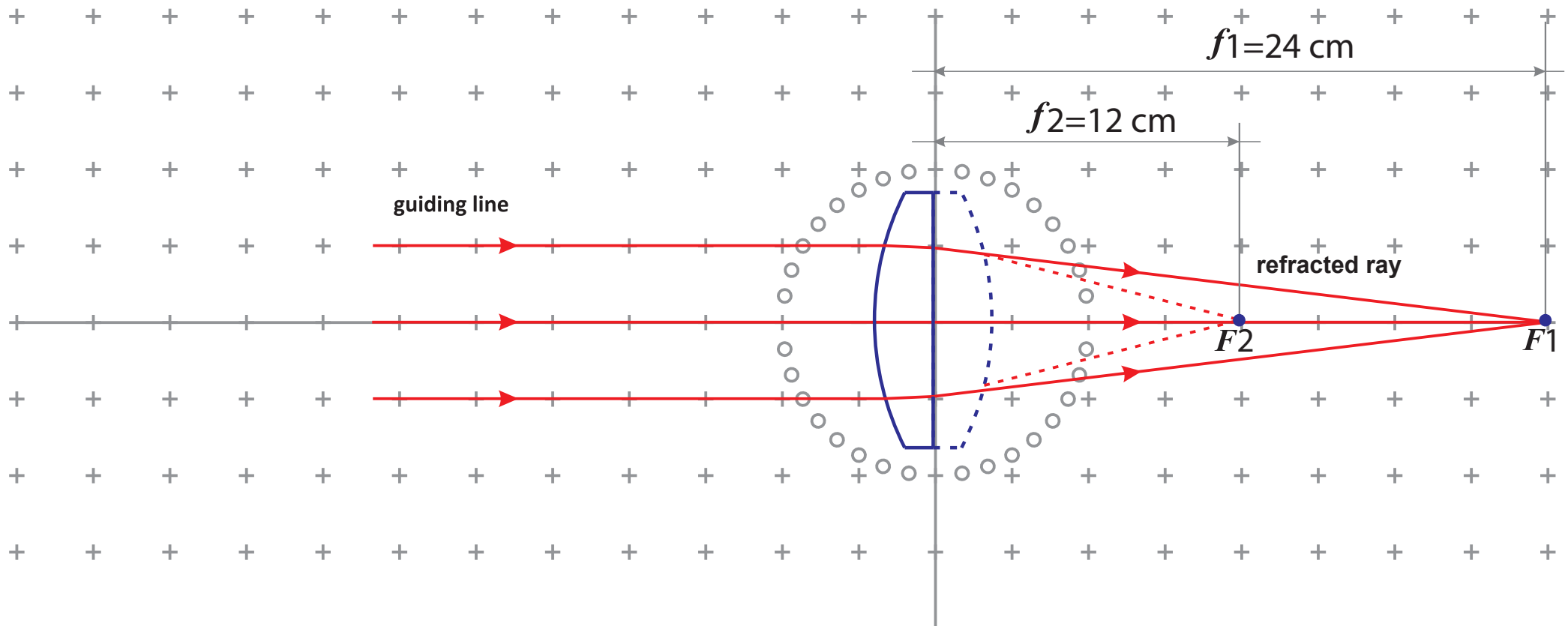

7- Convex Lens

4- Real Image (Magnification equal to unity)

- Place a convex lens on the determined position in the figure. Place two or three light sources on the board.
- Light rays pass the point **A1** and fall on the lens. We assume **A1** to be the **real source** of light rays.
- Light rays collimate at the point **A2** after refraction, and **A2** becomes a new light source.
- In this case, **A1** is called a **real source** and **A2** its **real image**. Why?

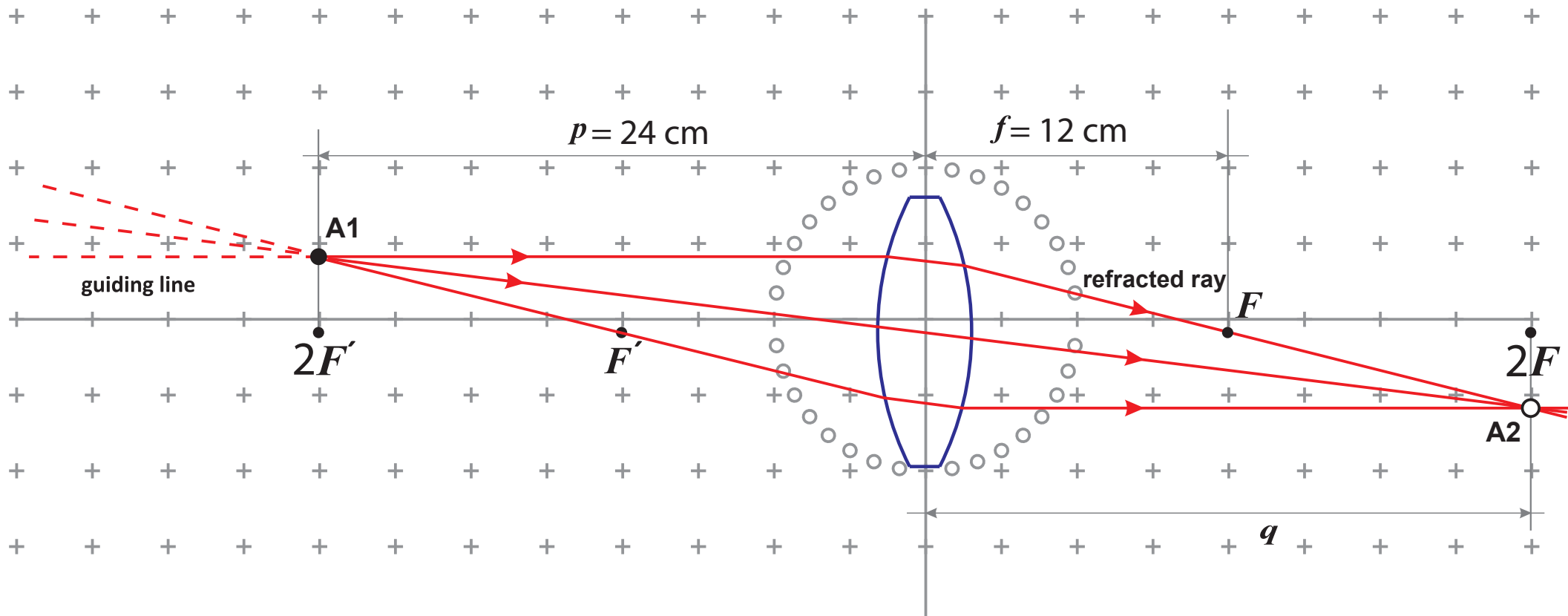


$$\frac{1}{p} + \frac{1}{q} = \frac{1}{f} \implies \frac{1}{24 \text{ cm}} + \frac{1}{q} = \frac{1}{12 \text{ cm}} \implies q = 24 \text{ cm}$$

7- Convex Lens

5- Real Image (Object: Infinite number of rays, Image: Infinite number of rays)

- Place the convex lens on the determined position in the figure. Place the mirrorlet on the point A1.
- Light ray falls on the mirrorlet. By oscillating the mirrorlet, emit light rays from A1 to the lens and get their refractions.
- A2 is the **real image** of A1. Repeat this experiment for other points of the optical object A1B1C1D1.
- Each **optical object and its image** are consisted of **infinite number of point sources** and from **each source** **infinite number of rays** is emitted.

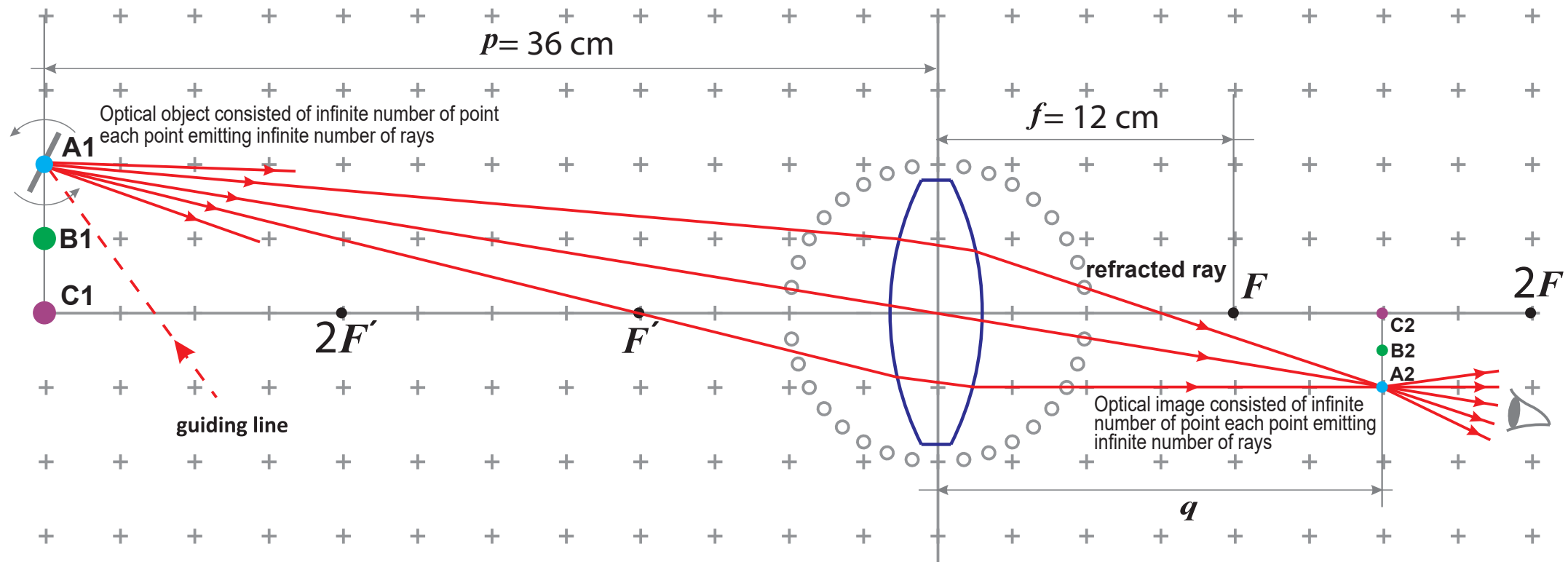


$$\frac{1}{p} + \frac{1}{q} = \frac{1}{f} \Rightarrow \frac{1}{36\text{ cm}} + \frac{1}{q} = \frac{1}{12\text{ cm}} \Rightarrow q = 18\text{ cm}$$

8- Concave Lens

1- Optical Center


- Place a concave lens on the determined position in the figure. Place two or three light sources on the board.
- With the aid of guiding lines set the incident rays such that they fall on the **optical center** of lens.
- Make sure that the lens position is precise, and that the light rays pass through the optical center.
- Each incident ray falling on the optical center, will pass through the lens **without refraction** and deviation in its path.

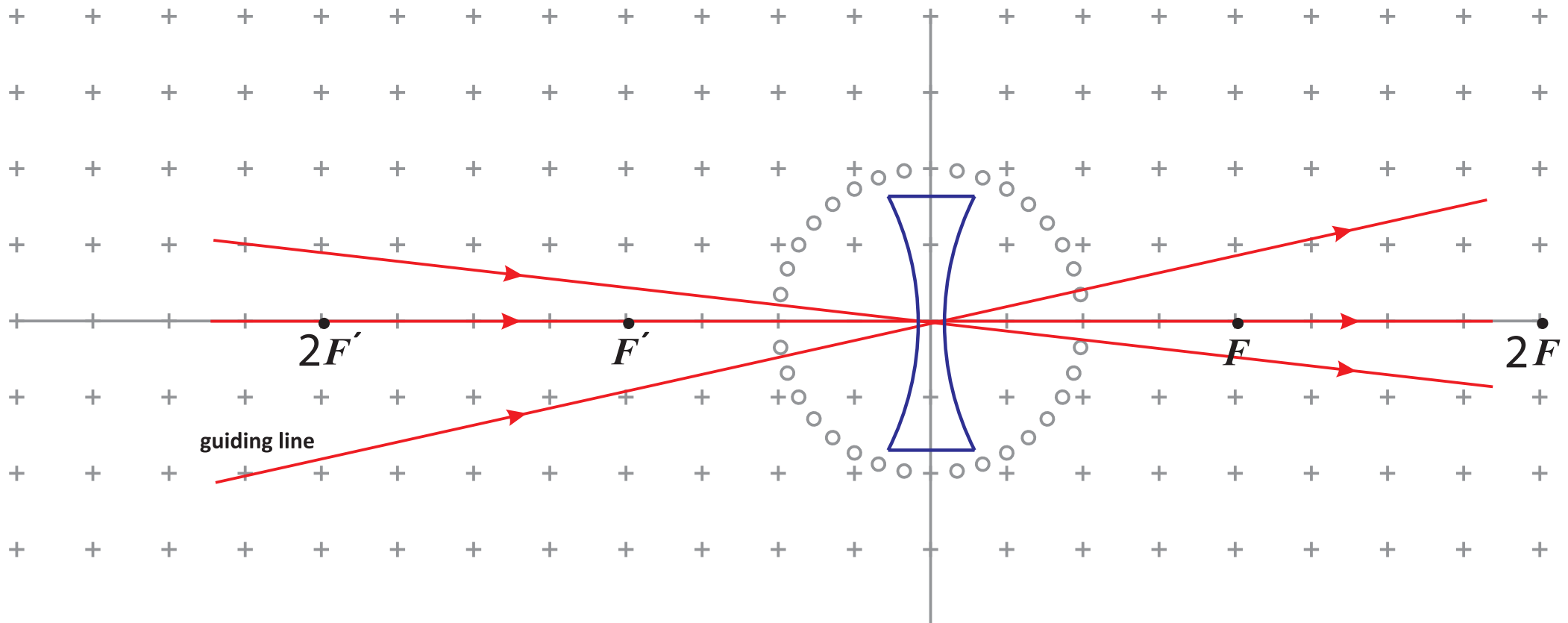

8- Concave Lens

2- Virtual Focal Point (Incident rays toward the virtual focal point become parallel)


- Place a concave lens on the determined position in the figure. Place two or three light sources on the board.
- With the aid of guiding lines emit light rays toward the focal point at the other side of concave lens.
- Make sure that the lens position is precise, and that the direction of incident rays pass the focal point.
- Incident ray toward the virtual focal point of the concave lens, will become parallel to the principal axis after refraction.

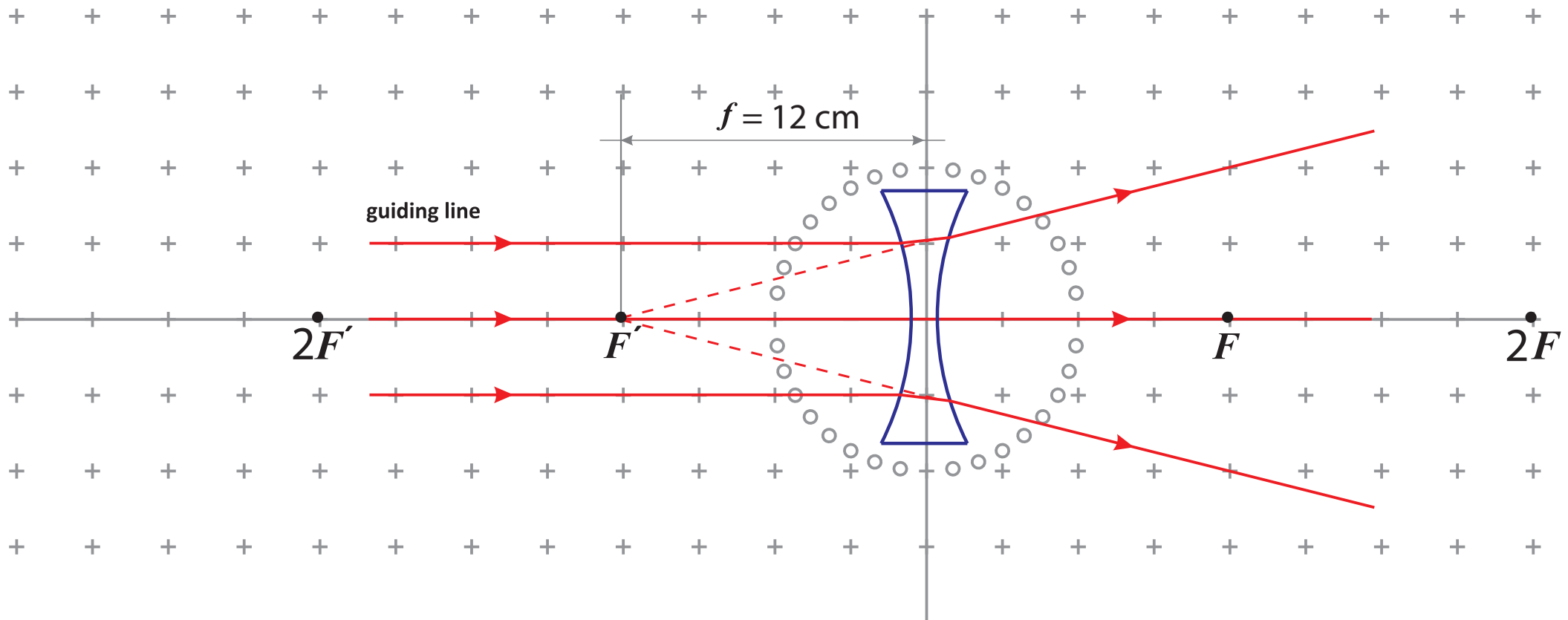

8- Concave Lens

3- Virtual Image (Magnification less than unity)

- Place a concave lens on the determined position in the figure. Place two or three light sources on the board.
- Light rays pass the point A1 and fall on the lens. We assume A1 to be the real source of light rays.
- Light rays after refraction diverge such that it seems they have emerged from the point A2.
- In this case A1 is called a **real source** and A2 its **virtual image**. Why?

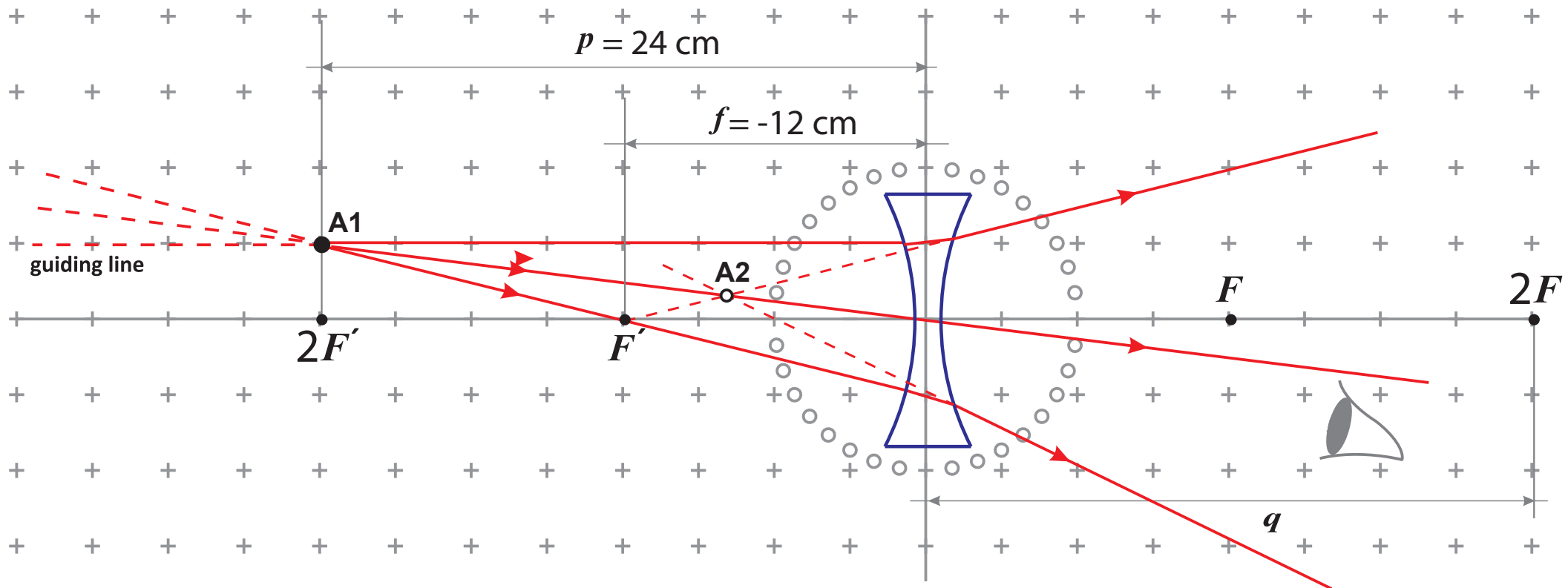


$$\frac{1}{p} + \frac{1}{q} = \frac{1}{f} \implies \frac{1}{24\text{ cm}} + \frac{1}{q} = \frac{1}{-12\text{ cm}} \implies q = -8\text{ cm}$$

9- Physics of the Eye

1- Healthy Lens

- The following figure is a schematic of the eye sphere. Place a convex lens suitable for the figure, as the **eye lens** on the determined position.
- By two or three light sources illuminate the lens with incident rays parallel to the principal axis.
- Make sure that the position of lens is precise and that light rays are parallel to the principal axis.
- Parallel rays are collimated on the eye's **retina**. It means that the image of a very distant point falls on the retina.

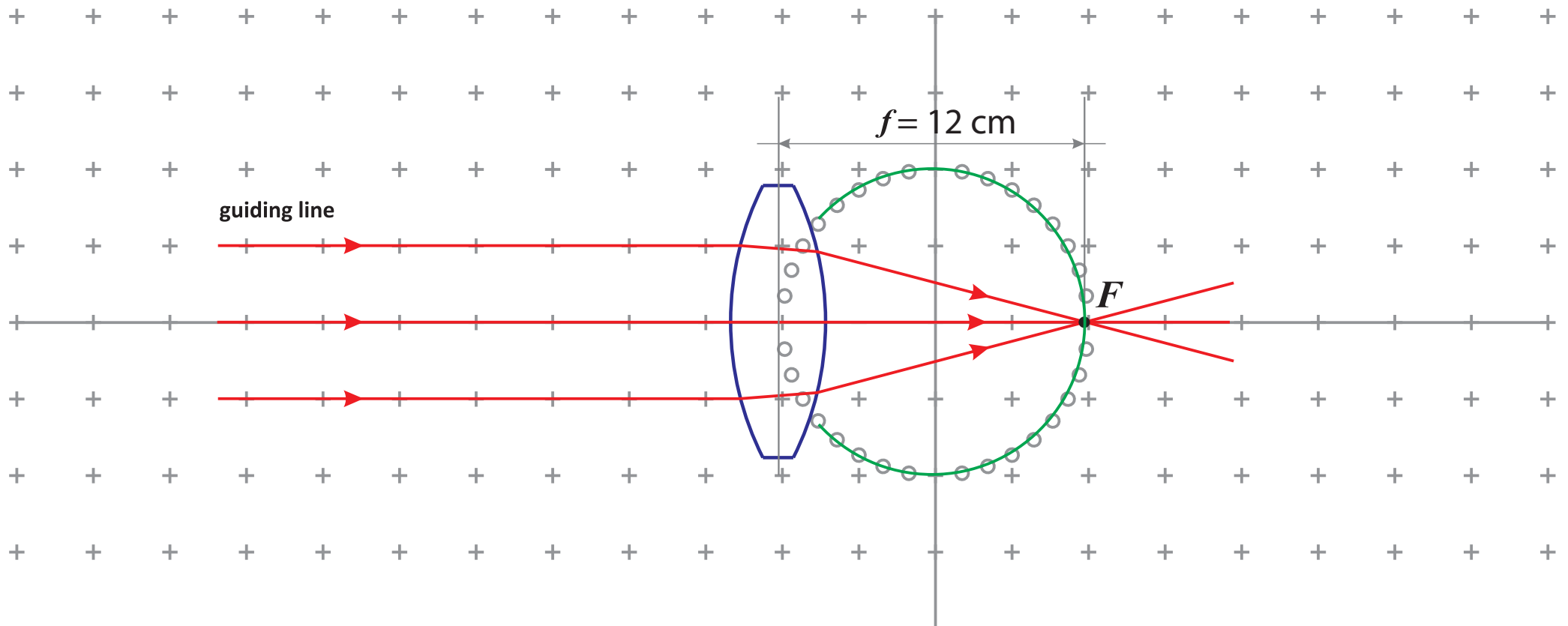

9- Physics of the Eye

2- Correction of the lens (Myopia (short-sightedness))

- The following figure is a schematic of the eye sphere. Place a convex lens suitable for the figure, as the **eye lens** on the determined position.
- By two or three light sources illuminate the lens with incident rays parallel to the principal axis.
- Rays are collimated in a position **in front of the retina**. This is a **short-sighted eye**.
- Solve the problem with a **concave lens** as an **eyeglass**.

